


BeadBug™-6 Quick Start Guide

Important! The BeadBug-6 is shipped with motor stabilizing screws in the base. These *must* be removed before use. Failure to do so may damage the unit.


1. Upon removing the unit from the carton, check to be sure that the voltage rating of the unit matches the local power supply and that the proper power cord has been received.
2. Locate the two shipping screws on the bottom of the unit (see image below). Using the supplied Allen key, remove these screws.

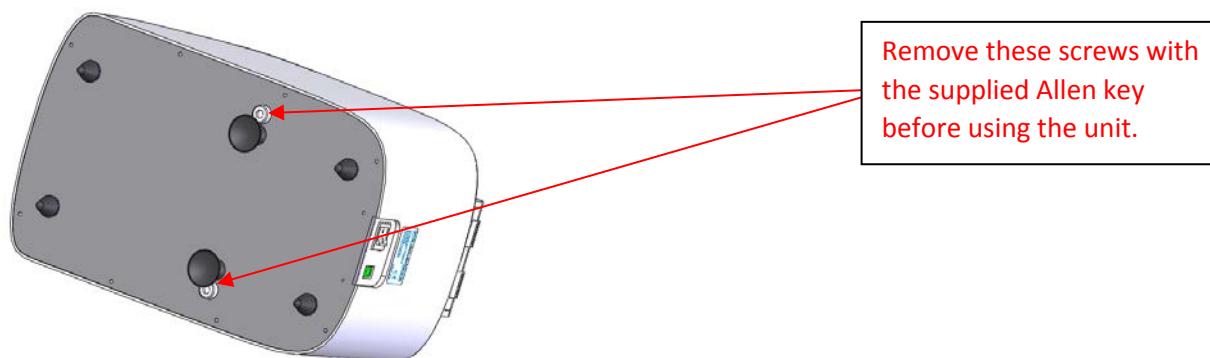

3. Connect the power cord to the outlet and switch on the power.
4. Load tubes symmetrically into the tube holder as shown in the illustration below. Do not run the unit without at least two tubes in place.



5. Press tubes firmly into place and lock the tube retaining bar into place as shown.

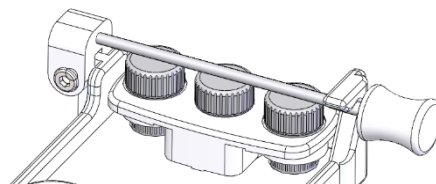

6. Be sure that the transparent lid is closed and locked in place.
7. Press "Set" followed by the up and down arrow keys to set the operating parameters. Pressing "set" additional times will scroll through the settable parameters: cycle number, run speed, run time and dwell time between cycles. Press "start" to begin the run
8. Should you need to stop the run, simply press "stop".

This Quick Start Guide is not intended to replace the operating manual. Please read the manual thoroughly before attempting to operate the BeadBug-6.