


MAGic Clamp™ Operating Instructions

To use the Magic Clamp, adjust the clamp to the appropriate flask size by following the instructions below:

- 1.) Press all three flask arms inward, to the inmost position
- 2.) Grip the sliding notch along the side wall of the clamp and turn until the pins are aligned with the inmost (125ml), middle (250ml) or outmost (500ml) position,
- 3.) Pull each of the three arms back until a “click” is heard, confirming proper positioning.

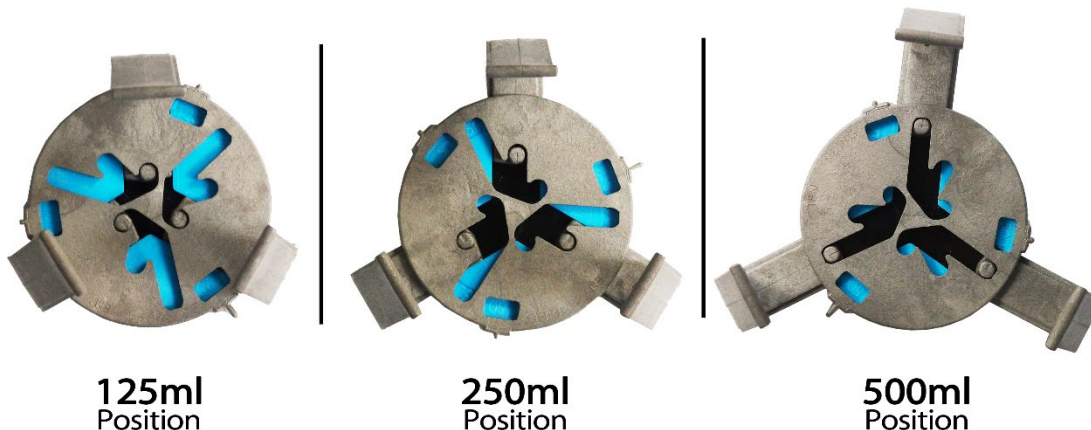

Once the clamp has been adjusted to the appropriate setting, place the clamp onto the **MAGic Clamp platform** in the desired platform location. The MAGic Clamp platform should consist of a steel base and a rubber mat. The platform is to be applied above the rubber mat.


Always attempt to align flasks in a balanced orientation. (Meaning, do not place all of the large sized flasks on one side of the platform and small sized flasks on the opposite side of the platform.)

MAGic Clamps are **ONLY** recommended for use with compatible MAGic Clamp platforms.

MAGic Clamp™ Operating Instructions

To use the Magic Clamp, adjust the clamp to the appropriate flask size by following the instructions below:

- 1.) Press all three flask arms inward, to the inmost position
- 2.) Grip the sliding notch along the side wall of the clamp and turn until the pins are aligned with the inmost (125ml), middle (250ml) or outmost (500ml) position,
- 3.) Pull each of the three arms back until a “click” is heard, confirming proper positioning.

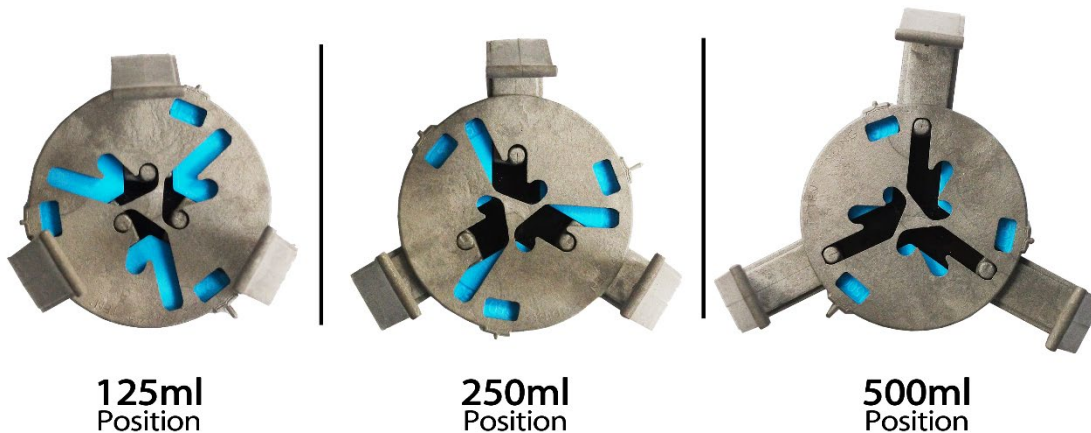

Once the clamp has been adjusted to the appropriate setting, place the clamp onto the **MAGic Clamp platform** in the desired platform location. The MAGic Clamp platform should consist of a steel base and a rubber mat. The platform is to be applied above the rubber mat.

Always attempt to align flasks in a balanced orientation. (Meaning, do not place all of the large sized flasks on one side of the platform and small sized flasks on the opposite side of the platform.)

MAGic Clamps are **ONLY** recommended for use with compatible MAGic Clamp platforms.