

## Setting up the Benchmark H2300-REG

1. Turn the knob on the CO2 tank to the closed position
2. Attach the large threaded adapter (B) to the nozzle of the tank with the white washer in between – FULLY TIGHTEN using an adjustable wrench (Tank nozzle should be a **CGA-320** size)
3. Attach the regulator (A) to the other end of the large threaded adapter – FULLY TIGHTEN using an adjustable wrench
4. Thread the desired hose adaptor (D or F) through the small threaded adapter (C or E) to the side output of the regulator – FULLY TIGHTEN using an adjustable wrench
5. Attach the black knob to the front of the regulator – only turn 1-2 times
6. Turn the knob on tank to the open position
7. SLOWLY turn the black knob on the front of the regulator clockwise to open. Slowly turn until the dial reaches 5 to 10 L/min.

