ST20D Pen DO meter Instruction Manual

Welcome to OHAUS (www.ohaus.com)!


After more than a century of perfecting the art of measurement through our durable weighing products, OHAUS precision is now available in a line of portable electrochemistry products that provides accurate measurements of DO, pH, ORP, conductivity, TDS and salinity.

ST series pen meters are the economical option when you are simply looking for meters that are easy to use and provide accurate measurements.

The pen meters should not be used by Children of age 12 or younger.

This product conforms to the RoHS Directive 2011/65/EC and the EMC Directive 2014/30/EU. The Declaration of Conformity is available online at www.ohaus.com/ce

Thank you for choosing OHAUS waterproof pen Dissolved Oxygen meter ST20D. Please read the manual completely before use.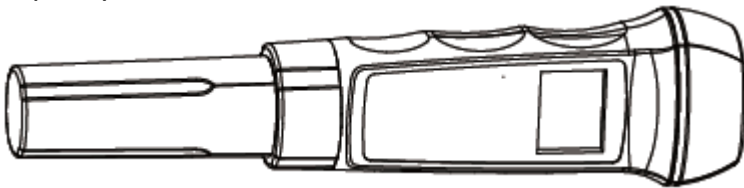

First Usage

Take off the membrane part, fill the electrolyte, then turn on the membrane. Warm up the DO pen meter by turning it on. The meter will turn off automatically after 10 minutes.

Calibration and measurement

After warm up, turn on the meter. Put the ST20D in the air and press the CAL button;

The meter will display "100" and blink under the signal, now the meter is calibrating 100% DO value; after 3 to 5 minutes when the signal is stable, press the Hold/Enter button to finish 100% DO calibration.

Now we have two options:

1. Press the Hold/Enter button to finish 1 point calibration and exit to measurement interface.

Note: Ohaus suggest that only do one point (100%) calibration for ST20D. Do not need perform 2 point (0%) calibration if the sample DO value is not too low (e.g. <1 mg/L).

2. Put the ST20D into the Zero-Oxygen solution. Press the CAL button to do the 0% point calibration, LCD will display and blink "0" under the signal, after about 3 minutes when the signal is stable, press the Hold/Enter button to finish 0% calibration.

Zero-Oxygen solution is saturated Na_2SO_3 (CAS NO 7757-83-7) solution. Ohaus offer DO zero-oxygen chemical (30059257).

After calibration, rinse the pen DO tip, then we can measure the sample.


Maintenance and Storage


After the usage, turn off the membrane, rinse the electrode and membrane, makes them dry and turn on the membrane part to store.

Harsh samples such as sticky, dirty or oily liquids may shorten the life time of the pen meter.


The meter uses 4 (1.5V AA) button batteries. Replace all batteries when the display fades or do not turn on.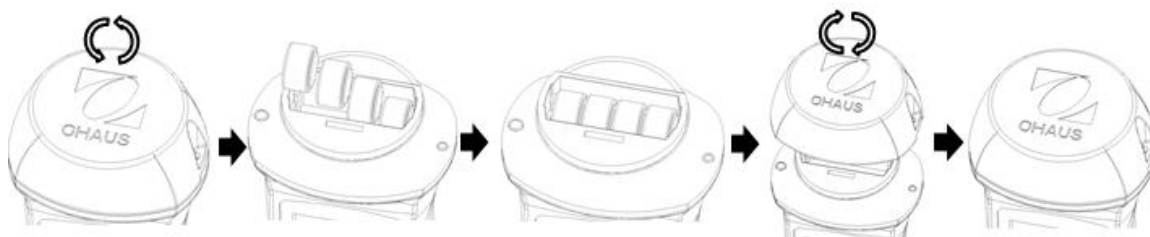

The electrode is replaceable.

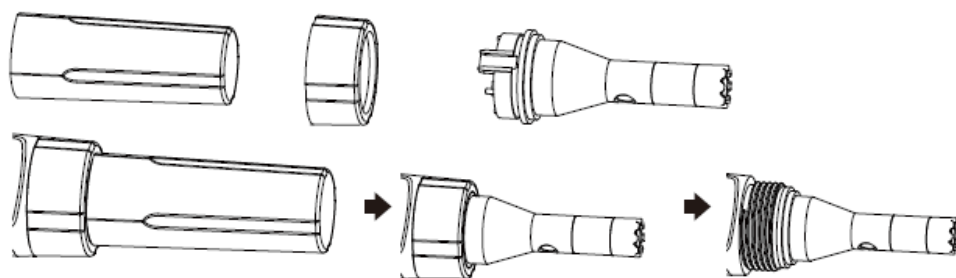The membrane part is replaceable.

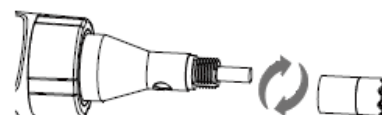Order Information

Model	Description	Item NO
ST20D	Polargraphic waterproof pen DO meter	30073986

Accessory

30222083	Pen Electrode, ST20D
30222084	Membrane part with electrolyte for ST20D
30059257	DO Zero Oxygen Chemical

Specification

Water Proof Pen DO Meter	ST20D
Measurement range	0.0 – 19.9 mg/L
Resolution	0.1 mg/L
Accuracy	± 0.5 mg/L
Battery	4 x 1.5V
Dimension	183mm x 45mm x 35mm
Weight	90g
Auto-off	After 10 min no operation

Auto temperature compensation range	5~45 °C (41-122 °F)
Materials	ABS

In conformance with the European Directive 2002/95/EC on Waste Electrical and Electronic Equipment (WEEE) this device may not be disposed of in domestic waste.

This also applies to countries outside the EU, per their specific requirements. Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. If you have any questions, please contact the responsible authority or the distributor from which you purchased this device.

Should this device be passed on to other parties (for private or professional use), the content of this regulation must also be related. For disposal instructions in Europe, refer to www.ohaus.com/weee. Thank you for your contribution to environmental protection.

FCC Notice: This equipment has been tested and found to comply with the limits for a Class B digital device, pursuant to part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference in a residential installation. This equipment generates uses and can radiate radio frequency energy and, if not installed and used in accordance with the instructions, may cause harmful interference to radio communications. However, there is no guarantee that interference will not occur in a particular installation. If this equipment does cause harmful interference to radio or television reception, which can be determined by turning the equipment off and on, the user is encouraged to try to correct the interference by one or more of the following measures: Reorient or relocate the receiving antenna, increase the separation between the equipment and receiver or consult the dealer or an experienced radio/TV technician for help.

The user is cautioned that any changes or modifications not expressly approved by the party responsible for compliance could void the user's authority to operate the equipment.

* 3 0 2 2 2 0 4 9 *

P/N 30222049 C ©2018 Ohaus Corporation, all rights reserved.

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