# PHCbi

## Operating Instructions

### Detector

# MLC-AD240A

**NOTICE:**

This product functions as a live cell metabolic analyzer in combination with the MLC-AC0 (controller). For proper handling and operation of the detector, please refer to controller's operating instructions in addition to these operating instructions.

Please read the operating instructions carefully before using this product and keep the operating instructions for future use.

See page 17 for the model number.

# CONTENTS

---

INTRODUCTION .....	3
SAFETY PRECAUTIONS.....	4
COMPONENTS .....	9
Detector components .....	9
Detector accessories.....	10
ROUTINE MAINTENANCE .....	11
Cleaning the detector .....	11
DISPOSAL OF UNIT .....	12
SPECIFICATION .....	17
SAFETY ENVIRONMENTAL CONDITIONS .....	17
SAFETY CHECK SHEET .....	18

# INTRODUCTION

---

- Read the operating instructions carefully before using the product and follow the instructions for safe operation.
- PHC Corporation takes no responsibility for safety if the product is not used as intended or is used with any procedures other than those given in the operating instructions.
- Keep the operating instructions in a suitable place so that they can be referred to as necessary.
- The operating instructions are subject to change without notice for improvement of performance or function.
- Contact our sales representative or agent if any page of the operating instructions is lost or the page order is incorrect, or if the instructions are unclear or inaccurate.
- No part of the operating instructions may be reproduced in any form without the express written permission of PHC Corporation.

## **IMPORTANT NOTICE**

PHC Corporation guarantees this product under certain warranty conditions. However, please note that PHC Corporation shall not be responsible for:

- any accident or damage caused by incorrect installment or handling of the detector

### <Intended Use>

This product is designed for the continuous measurement of the glucose concentration and lactate concentration in a culture medium during animal cell culture. This product cannot be used for diagnostic purposes.

# SAFETY PRECAUTIONS

Be sure to observe the operating instructions as they contain important safety advice. If the product is used in a manner not specified by the operating instructions, the protection provided by the product may be impaired.

For correct and safe use of the product, follow the precautions and procedures in these operating instructions carefully. Failure to do so could result in injury or damage to the product.

Precautions are illustrated in the following way:

## **WARNING**

Warning indicates a potentially hazardous situation which, if not avoided, could result in serious injury or death.

## **CAUTION**

Failure to observe CAUTION signs could result in injury to personnel and damage to the unit and associated property.

The following symbols are used in this document and some of them are attached to the unit.

	Actions are prohibited.
	Actions are mandatory.
	Caution must be taken.
	This symbol indicates a risk of an electric shock by touching the product with wet hands.
	This symbol indicates a risk of an electric shock by an electric leakage caused by a wet product.
	This symbol indicates a risk of injuries such as an electric shock caused by disassembling the product.
	This symbol indicates that the user must disconnect the mains plug for the purposes of maintenance, in the case of malfunction or when left unattended.
	This symbol indicates a potential danger or risk by biohazardous material. Ce symbole indique un danger ou un risque potentiel dû à des matières présentant un danger biologique.
	Read the operating instructions carefully before using the product.

---

**THE FOLLOWING APPLIES ONLY IN THE U.S.A.**

This device complies with part 15 of the FCC Rules. Operation is subject to the following two conditions:

- (1) This device may not cause harmful interference, and
- (2) This device must accept any interference received, including interference that may cause undesired operation.

**FCC Note:**

This product has been tested and found to comply with the limits for a Class A digital device, pursuant to part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference when the product is operated in a commercial environment. This product generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the instruction manual, may cause harmful interference to radio communications. Operation of this product in a residential area is likely to cause harmful interference in which case the user will be required to correct the interference at his own expense.

**THE FOLLOWING APPLIES ONLY IN CANADA.**

CAN ICES-003(A)/NMB-003(A)

# SAFETY PRECAUTIONS

## WARNING

### Installation

- **Do not install the detector at a place higher than your eye level.**  
The detector at a high place makes it difficult to locate the sensor module assembly on the tray, so that you may spill the culture medium over your body, causing an adverse effect on your body.

- **Before connecting the detector to the controller, turn off the controller and pull out the power plug from the outlet.**  
Otherwise, you may have an electric shock or get injured when you inadvertently touch an electric part.
- **Make sure that the detector cable and the connection cable are connected properly.**  
Otherwise, inadequate connection may cause a fire or an electric shock.
- **Handle the connection cables carefully.**  
Putting the controller or detector on them or bending/twisting them excessively may cause a short circuit or disconnection of a wire, which may lead to fire or an electric shock.
- **When you handle a harmful reagent, handle it in an isolated facility prepared for such operations.**  
Otherwise, an improper use may cause harm to the human body or natural environment.

### Connection cable

- **Do not handle the connection cable in a way that damages it (e.g., scratching, modifying, putting it close to a heat source, bending with force, twisting, pulling, adding weight, or binding).**  
Using the damaged connection cable may cause fire or an electric shock. Consult our services agency when you want to repair the connection cable.

- **Do not insert or pull out the connection cable with wet hands.**  
It may cause an electric shock.

- **Before cleaning, maintenance, or checkup, disconnect the connection cable.**  
Otherwise, you may have an electric shock or injury.
- **When disconnecting the connection cable, hold the connector at the end of the cable.**  
Pulling the cable may cause an electric shock.

# **WARNING**

## **When using the product**

- **Protect from strong vibration or impact when moving or operating the detector.**  
The detector may be damaged and cause a fire.
- **Do not put any foreign objects such as a pin or metal piece into any gap or hole of the detector, or the inside of the detector when the detector tray is pulled out.**  
It may cause an electric shock or injury by accidental contact with moving parts.

- **Do not use the detector when a foreign object or liquids such as reagents have spilled into the detector or when the detector fell to the ground.**  
If you continue to use it under a defective condition, it may cause a fire or other problems. Contact our service office or sales agency right away for inspection and repair.

- **Do not spill water directly over the detector.**  
Spilled water may cause an electric shock or fire.

- **When you open or close the detector tray, open or close it carefully.**  
If you open or close the tray quickly, the culture medium may overflow inside the sensor module assembly and enter inside the detector, causing contamination.
- **When installing the detector in the CO<sub>2</sub> incubator, do not put any objects on trays set above the detector.**  
If a culture medium spills, it may enter inside the detector, causing contamination or a fault.

## **When using a reagent and samples**

- **Wear appropriate protective gear such as gloves and glasses when you handle hazardous substances such as lactate solution.**  
A lactate solution causes a serious injury if it touches your skin or eye.

- **Observe the laboratory biosafety guidelines issued by WHO.**  
In addition to the precaution described here, refer to the laboratory biosafety guidelines issued by WHO. This product is assumed to be used in a laboratory at biosafety level 2 or lower.
- **Wear appropriate protective gear when handling a potentially infective sample or a product that may had contact with such a sample.**  
Touching them directly may cause infection.
- **Clean the parts that might have come into contact with a potentially infective sample.**  
An infective sample left on the surface of the parts may cause infection.

# SAFETY PRECAUTIONS

## WARNING

### When something is wrong with the product

- **When an error occurs, immediately turn off the controller and pull out the power plug.**  
If you continue to use it under a defective condition, it may cause a fire or other problems. Contact our service office or sales agency right away for inspection and repair.

- **Do not disassemble or modify the detector.**  
It may cause an electric shock. Contact our service office or sales agency for the inspection or repair.

### When storing and disposing of the product

- **If the product will not be used for a long time, remove the detector from the CO<sub>2</sub> incubator, and store it.**  
Otherwise, the deterioration of insulation may cause an electric shock, electric leakage, or fire.

- **Ask a qualified contractor to disassemble or dispose of the products. Do not leave the unit in a location that can be accessed by third parties.**  
Otherwise, an unexpected accident (by unintended use, etc.) may happen.

- **Observe the ordinances or regulations of the local government when disposing of the products that had contact with an infectious sample.**  
Otherwise, it may cause infection.

## CAUTION

- **Be sure to wear gloves when you clean the detector.**  
Otherwise, you may be injured by the parts inside the detector.
- **Do not insert your hand in the opening of the detector tray.**  
You may be injured by the parts inside the detector.
- **When you carry the detector, carry it with both hands.**  
Otherwise, you may be injured when you drop it.
- **Confirm that the CO<sub>2</sub> incubator tray is strong enough to bear the detector.**  
Otherwise, the tray may collapse to cause injury.

- **Before moving the detector, take the sensor module assembly out of the detector.**  
Otherwise, the sample may spill within the detector, contaminating the detector. Also, a spilled sample may pass infection to humans.
- **Make sure to clean the detector before storing or transporting it.**  
A contaminated detector may cause infection to humans.


- **Be careful not to pinch your fingers when you close the detector tray.**  
Otherwise, you may be injured.

# COMPONENTS

## Detector components

### Front and tray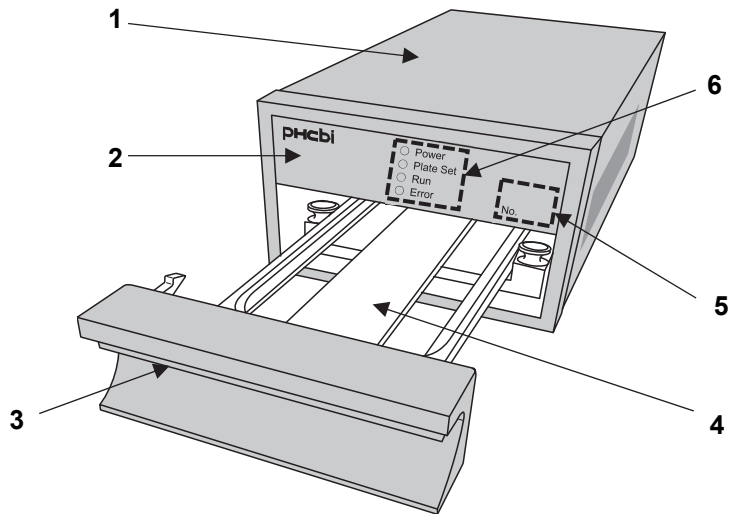

#### 1. Housing

The housing covers the detector. There are ventilation openings on the sides of the housing. Do not block the ventilation openings when you install the detector.

#### 2. Front panel


The detector status and detector ID number are indicated on the front panel.

#### 3. Tray trigger

Pull the tray trigger toward you to draw the tray out.

#### 4. Tray

Put the sensor module assembly (or check module assembly) on the tray.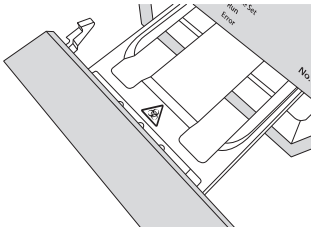

 <b>WARNING</b>	<p>The sensor module assembly may include a potentially infective sample. Clean the tray when such a sample has been attached to the tray.</p> <p>An infective sample left on the tray may cause infection.</p>
--	---

#### 5. Detector ID number indication

Displays the detector ID number that is recognized by the controller.

#### 6. Status LED

Indicate the detector status.


Status LED	Condition	Details
Power (green)	Lighting	The detector is connected to the controller that is turned on.
Plate Set (white)	Lighting	The sensor module assembly (or check module assembly) is set on the tray in the detector.
Run (white)	Blinking	A voltage is being applied to the sensor module.
	Blinking (fast)	The assay is paused.
	Blinking (fast)	A plate set error has occurred.
Error (red)	Blinking (fast)	An error that needs to be recovered by user operation has occurred.
	Lighting	An error that does not need to be recovered by user operation has occurred.

#### Note:

After the controller is turned on, all status LEDs are lit for five seconds when the detector is recognized by the controller.

# COMPONENTS

## Side, rear, and bottom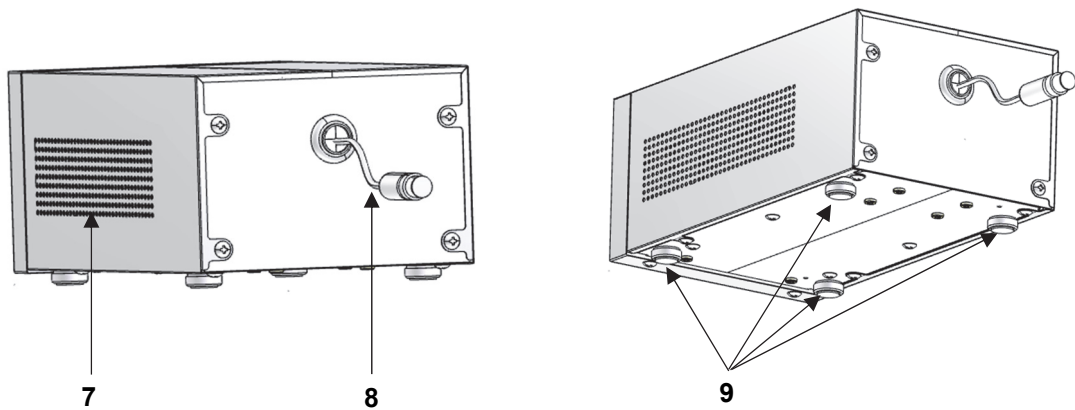

**7. Ventilation openings (on the left and right sides)**


Do not block them when you install the detector. If the ventilation openings are blocked, CO<sub>2</sub> gas and humidity do not circulate and may affect cultivation.

**8. Detector cable**

Connect it to the connection cable for connecting the detector and the controller. The connector of this cable is waterproof.



**9. Leveling feet**

Turn the leveling feet to adjust the lengths of legs. Use the four leveling feet to set the detector horizontally.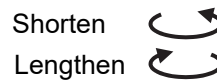

## Detector accessories

Check that the following accessories are included with the detector. If anything is missing, contact our sales representative or agent.

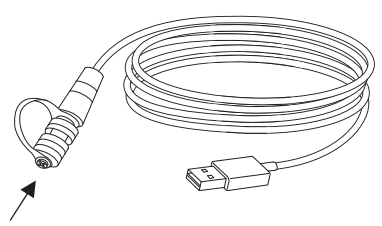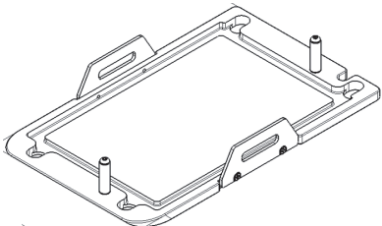Name	Quantity	Appearance
Connection cable (with a protection cap)	1	 <p>Be sure to cover the waterproof connector with the protection cap when the detector is not connected.</p>
Plate adapter (bottom)	1	

# ROUTINE MAINTENANCE

In order to obtain precise data, it is necessary to clean the products regularly. Follow the sections below to clean the detector. Also, yearly maintenance by the manufacturer is recommended. Contact our sales representative or agent for maintenance requests.


## Notes:

- Clean the exterior of the detector, detector tray, and cables using 70% ethanol once a month.
- Contact our sales representative or agent to disassemble and clean the inside of the detector once a year. Do not take apart the detector yourself because it may lead to an accident or failure of the detector.
- Contact our sales representative or agent when the products are too dirty.
- Do not use brushes, acid, alkaline or chlorine-based detergents, disinfectants, volatile chemicals such as thinner, soap powder, polishing powder (cleansers), or boiling water. It may cause discoloration, corrosion, and rust.
- Do not clean the products unit using peracetic acid or hydrogen peroxide, or do not disinfect the products by formalin fumigation.
- Never pour water onto or into the products. Doing so can damage the electric insulation and cause failure.
- Turn off the product before cleaning.

## Cleaning the detector

**CAUTION** Do not put your hands into the inside of the detector when the tray is pulled out. You may get injured by touching the interior parts.

1. Turn off the power to the CO<sub>2</sub> incubator and take the humidifying pan out of the incubator by following the operating instructions for the CO<sub>2</sub> incubator.
2. Put on rubber gloves, and then disinfect the surface of the rubber gloves with 70% ethanol.
3. Open the incubator doors to cool down the incubator and reduce humidity in the chamber. Wipe the housing, front panel, tray trigger, and detector cable using a gauze moistened with a proper amount (the amount that cannot form droplets) of 70% ethanol. Do not spray 70% ethanol on the detector directly.
4. Fully pull out the tray of the detector and thoroughly wipe the area shown in the following figure.

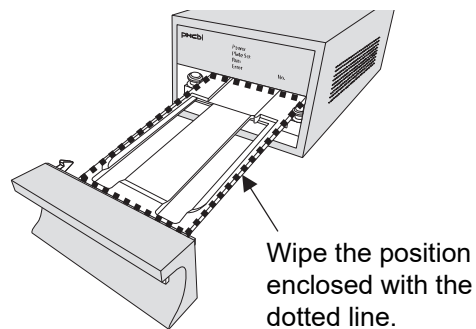# STORING THE PRODUCT

---

When the product is not used for a long time, store it by following the steps below. Storing the detector for a long time in the experimental condition may cause failure.

Before storing the detector, take the detector out of the incubator and clean it.

**Note:**

Check that the sensor module assembly (or check module assembly) is not left inside the detector. If the sensor module assembly is left in the detector, take it out of the detector.

	<b>Caution</b>	<b>Wear gloves and a mask during work.</b> Without them, you may be injured by edges of internal parts. Contacting chemicals and inhaling dust are bad for your health.
---	----------------	---

1. Wear rubber gloves and sterilize their surfaces with 70% ethanol.
2. Confirm that the temperature and humidity in the CO<sub>2</sub> incubator have dropped and turn off the incubator. Then, take the humidification pan out of the incubator and dispose of the water in the pan.
3. Pull out the power plug for the controller AC power cable from the outlet (turn off the detector).
4. Pull the waterproof connector connecting the detector and the controller to the front side.
5. Turn the outer ring of the waterproof connector counterclockwise to remove the connector.
6. Hold the detector with both hands tightly and take it out of the incubator.
7. Clean the detector by following the steps for cleaning (page 11).
8. After the ethanol has evaporated completely, put the detector in the plastic bag that came with the detector and store the detector (if you cannot find the plastic bag, use a clean store-bought plastic bag for storage).

# DISPOSAL OF UNIT

---

When disposing of the products, ask a qualified contractor.

	<b>WARNING</b>	<b>Ask a qualified contractor to carry out disassembly/disposal of the products and do not leave the product in a location that can be accessed by third parties.</b> This may result in unexpected accidents (e.g., the products may be used for unintended purposes).
---	----------------	--

**Notes:**

- Before disposing of the products, decontaminate the products to the extent possible by the user.
- For details on risks during storage, transportation, and disposal of the products, contact our agent listed on the separate information list.

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# SPECIFICATION

---

Product name	Detector
Model number	MLC-AD240A-PW
Dimension	Width: 162 mm, Depth: 290 mm, Height: 118 mm
Weight	4.7 kg
LED display	Status LED: POWER (green), PlateSet (white), RUN (white), and ERROR (red) Detector ID number indication: For displaying detector ID number
Applied voltage	RE voltage: 700 mV, WE voltage: 800 mV
Power source	USB bus power: 5 V DC
Maximum current consumption	65 mA
Environmental condition	Temperature: 37°C, CO <sub>2</sub> concentration: 5%, Humidity: 95% RH ± 5% RH *Expected to be used in a CO <sub>2</sub> incubator. Operation in other environments is not guaranteed.
Recommended number of units to be installed in an incubator	MCO-50: Max. 1 unit MCO-170: Max. 4 units* MCO-230: Max. 4 units* *When placing two detectors on one shelf, use a reinforced shelf. <b>Note:</b> When using a CO <sub>2</sub> incubator other than ours, the performance of the detector and incubator may be affected. If you wish to install the detectors in an incubator other than ours, please check the impact on the performance by yourself in advance.
Cleaning	Wipe with a cloth moistened with 70% ethanol. Hydrogen peroxide sterilization and dry-heat sterilization are not supported.
Accessories	Connection cable, plate adapter (bottom)

**Note:**

The product data has been measured based on the PHC standard.

# SAFETY ENVIRONMENTAL CONDITIONS

---

This equipment is designed to be safe at least under the following conditions (based on the IEC 61010-1):

- Indoor use;
- Altitude up to 2,000 m;
- Applicable pollution degree of the intended environment (POLLUTION DEGREE 2);

# SAFETY CHECK SHEET

**CAUTION**

Please copy and fill out this form before servicing.  
Hand over the form to the service engineer for their  
and your safety.

## Safety check sheet

### 1. Stored material

Risk of infection: Yes No Maybe

Risk of toxicity: Yes No Maybe

Risk from radioactive sources: Yes No Maybe

List all potentially hazardous materials that have been stored in this unit:

---

---

### 2. Contamination in the unit

a) Contamination Yes No Maybe

Types of contamination (if any): \_\_\_\_\_

---

b) Decontaminated Yes No

Methods used for the decontamination work:

---

---

### 3. Status of the unit

a) The unit is now safe to work on Yes No

b) If the answer is "No,"

Details on the danger: \_\_\_\_\_

---

Measures we should take to reduce the danger:

---

---

Date:

Signature:

Address, Division:

Telephone:

Product name:

Detector

Model No.

MLC-AD240A

Serial number:

Date of Installation:

Please decontaminate the unit yourself before calling the service engineer.

# MEMO

---

Manufacturer's name and address

**PHC Corporation**

1-1-1 Sakada, Oizumi-machi, Ora-gun, Gunma 370-0596, Japan

Importer's name and address

within the USA : **PHC Corporation of North America**

1300 Michael Drive, Suite A, Wood Dale, IL 60191

Toll Free USA (800) 858-8442

Importer's name and address

within the European Union : **PHC Europe B.V.**

Eikdonk 1, 4825 AZ, Breda, The Netherlands

within the United Kingdom : **PHC Europe B.V.**

9 The Office Village, North Road, Loughborough LE11 1QJ, United Kingdom

© PHC Corporation 2024

LDCL084400-0

Printed in Japan

LDCL084400-0

J0524-0