



Operating Instructions

Pharmaceutical Refrigerator

MPR-S1201XH-PA MPR-S1201RXH-PA



MPR-S1201XH-PA

Please read the operating instructions carefully before using this product, and keep the operating instructions for future use.

See page 63 for model number.

CONTENTS

INTRODUCTION	4
PRECAUTIONS FOR SAFE OPERATION	5
ENVIRONMENTAL CONDITIONS	10
SYMBOLS ON UNIT	11
REFRIGERATOR COMPONENTS	
Main body	12
Control panel	14
INSTALLATION SITE	16
INSTALLATION	17
PLACEMENT OF STORED ITEMS	19
USE OF SHELVES AND DRAWERS	20
START-UP PROCEDURE	21
OPERATION DURING POWER FAILURE	21
AFTER RECOVERY FROM POWER FAILURE	21
HOME SCREEN (Temperature display screen)	22
MENU SCREEN	23
SETTING THE CHAMBER TEMPERATURE	25
SETTING THE KEYPAD LOCK	26
RELEASING THE KEYPAD LOCK	27
DISPLAY OF MAX/MIN TEMPERATURE	28
RESET OF MAX/MIN TEMPERATURE	29
SETTING THE PERIOD FOR MAX/MIN TEMPERATURE	30
DISPLAY AND EXPORT OF DATA LOG	
Display & export of temp. and door operation logs	31
Export of temperature and door operation logs	33
Setting the log interval	34
Setting the device ID	35
DISPLAY OF EVENT LOG	36
EXPORT OF EVENT LOG	37
SETTING THE SELECTABLE PARAMETERS	
Selectable parameters	38
Setting the temperature display	39
Setting the high-temperature alarm	40
Setting the low-temperature alarm	41
Setting the door buzzer delay	42
Setting the buzzer suspended period	43
Setting the buzzer volume	44
Setting the remote alarm	45
Setting the interior light	46
Setting the lighting time	47
Setting the compressor delay	48

CONTENTS

Setting the rail heater	49
Setting the date display (format)	50
Setting the date	51
Setting the time	52
REMOTE ALARM TERMINAL	53
ALARMS AND SELF-DIAGNOSIS	54
ROUTINE MAINTENANCE	
Cleaning of exterior, interior, door rail (including drain port) and accessories	56
DEFROSTING	57
TROUBLESHOOTING	58
DISPOSAL OF UNIT	60
Decontamination of unit	60
TEMPERATURE RECORDER (OPTION)	61
BATTERY KIT FOR POWER FAILURE ALARM (OPTION)	61
SPECIFICATIONS	62
PERFORMANCE	63
SAFETY CHECK SHEET	64

INTRODUCTION

Read the operating instructions carefully before using the appliance and follow the instructions for safe operation.

Handling of Operating Instructions

- PHC Corporation never guarantees any safety if the appliance is used for any objects other than intended use or used by any procedures other than those mentioned in the operating instructions.
 - Keep the operating instructions in a suitable place to refer to it as necessary.
 - Contact our sales representative or agent if any page of the operating instructions is lost or page order is incorrect.
 - Contact our sales representative or agent if any point in the operating instructions is unclear or if there are any inaccuracies.
- ◇The operating instructions is only for the pharmaceutical refrigerator MPR-S1201XH and MPR-S1201RXH manufactured by PHC Corporation.
- ◇The contents are subject to change without notice.
- ◇No part of the operating instructions may be reproduced in any form without the expressed written permission of PHC Corporation.

Proposal for safe storage of refrigerator contents

We can offer some precautions below to protect the valuable material against accidental temperature rise. Contact our sales representative or agent for the details or installation.

- Temperature recorder (Option)
- Battery kit for power failure alarm (Option)
- Remote alarm system (Marketed product)

IMPORTANT NOTICE

PHC Corporation guarantees this product under certain warranty conditions. However, please note that PHC Corporation shall not be responsible for any loss or damage to the contents of the product.

<Intended Use>

This equipment is designed for storage of pharmaceuticals, samples and reagents.

PRECAUTIONS FOR SAFE OPERATION

It is imperative that the user complies with the operating instructions as they contain important safety advice.

Items and procedures are described so that you can use this unit correctly and safely. Following these precautions will prevent possible injury to the user and any other person.

Precautions are illustrated in the following way:




WARNING

Warning indicates a potentially hazardous situation which, if not avoided, could result in serious injury or death.

CAUTION

Failure to observe CAUTION signs could result in injury to personnel and damage to the unit and associated property.

Symbols have the following meanings:

-  This symbol means caution.
-  This symbol means an action is prohibited.
-  This symbol means an instruction must be followed.











Be sure to keep the operating instructions in a place that is accessible to users of this unit.

For the State of California, USA Only:

This product contains a CR Coin Cell Lithium Battery which contains Perchlorate Material – special handling may apply. See www.dtsc.ca.gov/hazardouswaste/perchlorate.














PRECAUTIONS FOR SAFE OPERATION

WARNING

-  **Never damage the power supply cord or power supply plug (by breaking, adapting, placing near a source of heat, bending with force, twisting, pulling, adding weight, or binding).** A damaged power supply cord or power supply plug may cause electric shock, short circuit, or fire.
-  **Remove dust from the power supply plug periodically.** Dust on the power supply plug may cause insulation failure due to moisture and thus cause a fire. Disconnect the power supply plug and wipe it with a dry cloth.
-  **Make sure the power supply plug is pushed fully in.** Faulty insertion of the power supply plug may cause electric shock or fire due to generation of heat. Never use a damaged power supply plug or loose power outlet.
-  **Do not touch any electrical parts (such as power supply plug) or operate switches with a wet hand.** This may cause electric shock.
-  **Use a dedicated power source as indicated on the rating label attached to the unit.** Multiple socket adaptor may cause fire resulting from abnormal heating.
-  **Do not use the unit outdoors.** Exposure to rain may cause leakage and/or electric shock.
-  **Only qualified engineers or service personnel should install the unit.** Installation by unqualified personnel may cause water leakage, electric shock, or fire.
-  **Install the unit in a location capable of bearing the total combined weight (product + optional accessories + stored items). After installing the unit, be absolutely sure to take precautions to prevent the unit from falling over.** If the unit is installed in a location which is not strong enough or if the proper precautions are not taken, the unit may fall over and cause injuries.
-  **Do not install the unit in a location where flammable or volatile substances are present.** Installing the unit in a location where flammable or volatile substances are present may cause explosions and/or a fire.
-  **Do not install the unit where there are high levels of moisture or where it may be splashed with water.** Installing the unit where there are high levels of moisture or where it may be splashed with water may cause the insulation to deteriorate and give rise to leakage and/or electric shock.
-  **Be absolutely sure to earth (ground) the unit in order to prevent electric shock.** Failure to earth the product may give rise to electric shock. If necessary, ask a qualified contractor to do this work.
-  **Do not connect the earth wire to a gas pipe, water pipe, or lightning rod when earthing the unit.** Earthing the unit improperly may give rise to electric shock.












PRECAUTIONS FOR SAFE OPERATION

WARNING

-  **Never store volatile or flammable substances in this unit except in a sealed container.** Such substances may cause explosion or fire if they leak.
-  **Never insert metal objects such as pins and wires into any vent, gap, or outlet on the unit.** This may cause electric shock or injury by accidental contact with moving parts.
-  **When handling harmful samples (for example, those which consist of toxic, pathogenic, or radioactive substances), install the unit inside a designated isolation facility.** If the unit is installed in a location which is not an isolation facility, there may be detrimental effects on both people and the natural environment.
-  **Before proceeding with maintenance or checking of the unit, disconnect the power supply plug.** Performing the work while power is still flowing to the product or while the power supply plug is still connected may give rise to electric shock and/or injury.
-  **Wear protective gloves and mask during maintenance.** Touching or inhaling chemicals or aerosols from around the unit may be detrimental to health.
-  **Never splash water directly onto the unit** as this may cause electric shock or short circuit.
-  **Never put containers with liquid on top of the unit** as this may cause electric shock or short circuit if the liquid is spilled.
-  **Never disassemble, repair, or modify the unit yourself.** A high-voltage area is located inside the unit. Any work carried out by unauthorized personnel may result in electric shock. Contact our sales representative or agent for maintenance or repair.
-  **Disconnect the power supply plug if there is anything wrong with the unit.** Continued abnormal operation may cause electric shock or fire.
-  **Grip the power supply plug when disconnecting the power supply cord from the outlet.** Pulling the power supply cord may cause electric shock or short circuit.
-  **Disconnect the power supply cord when the unit is not in use for long periods.** Keeping the unit connected may cause electric shock, leakage, or fire due to the deterioration of insulation.
-  **Do not place this unit in a location where it is difficult to disconnect the power supply plug.** Failure to disconnect the power supply plug may cause fire in the event of a problem or malfunction.
-  **If the unit is to be stored unused in an unsupervised area for a long period, ensure that children do not have access and that doors cannot be closed completely.**










PRECAUTIONS FOR SAFE OPERATION

WARNING

-  **Ask a qualified contractor to carry out disassembly and disposal of the unit.** Leaving the unit in a location that can be accessed by third parties may result in unexpected accidents (e.g. the unit may be used for unintended purposes).
-  **Do not leave the plastic bags used for packing in a place where they can be reached by small children** as this may result in unexpected accidents such as suffocation.
-  **When moving the unit, be sure to take precautions to prevent it from falling over.** Moving the unit with too much force may cause it to fall over, possibly resulting in injury. A qualified individual must be assigned to supervise the safe movement and relocation of the unit.
-  **Do not install the unit in a location where corrosive gases such as acids are present.** Installing the unit in a location where corrosive substances are present may cause electric components to corrode, leading to leakage and/or electric shock due to the deterioration of insulation resulting from corroded electrical components.
-  **Connect the unit to a power source as indicated on the rating label attached to the unit.** Use of any other voltage or frequency other than that on the rating label may cause fire or electric shock.
-  **Never damage the chamber wall or pipework in the chamber when removing frost.** The refrigerant is flammable and may cause a fire if it leaks.
-  **Install the unit in a well-ventilated (airy) location to prevent the accumulation of flammable refrigerant.** The flammable refrigerant may cause fire if it leaks.
-  **Flammable and explosive product.** The unit contains flammable refrigerant. When repairing or recycling, only trained service personnel will repair and follow the procedure below.
 - Ensure the room is well ventilated to prevent refrigerant accumulation.
 - Keep fire away when the refrigerant is contained in the product.
 - Do not damage or break the pipework.
-  **Stop using the fire, open the windows for ventilation, and leave the unit as it is if the pipework is damaged or the leakage of flammable refrigerant is noticed. Then contact our sales representative or agent.** Disconnection or connection of the power supply plug may cause ignition, explosion, or fire.
-  **Always use the power supply cord that is provided.** Other power supply cord may cause electric shock or fire.
-  **Do not use the provided power supply cord for other equipment.** That may cause electric shock or fire.

PRECAUTIONS FOR SAFE OPERATION

CAUTION

-  **This unit must be plugged into a dedicated circuit** protected by branch circuit breaker.
-  **Use a dedicated power source as indicated on the rating label attached to the unit.** Multiple socket adaptor may cause fire resulting from abnormal heating.
-  **Do not climb on top of the unit or put any objects on the unit.** Falling from the unit may cause injury; falling objects may cause damage to the unit.
-  **Never install the unit in a location where corrosive materials such as sulphur compounds are likely to be generated (e.g. near a drainage facility).** Corrosion of the copper pipes may result in the deterioration and consequently the failure of the cooling unit.
-  **Before resuming operation after the unit has been turned off, be absolutely sure to check the settings.** Items stored inside the unit may be adversely affected if operation of the unit is resumed with changed settings.
-  **Never store corrosive substances such as acids or alkalis in the unit except in a sealed container.** These may be detrimental to health and may cause corrosion of internal components, cooling circuit or electrical components.
-  **To ensure the safety of the service engineer, submit a safety check sheet with the required items filled out.** This is provided as the photocopyable “Safety Check Sheet” at the end of these operating instructions.
-  **Use designated parts for parts replacement.** Using an incorrect part may cause fire.
-  **Do not cause strong shock or vibration during movement or use.** The piping may be damaged, causing a fire.
-  **Flammable and explosive product.** The unit contains flammable refrigerant. Consult repair manual/owner’s guide before attempting to install or service this product. All safety precautions must be followed.
-  **Flammable refrigerant used. Flammable and explosive product.** Dispose of properly in accordance with national regulations.
-  **Take care when handling the glass door when removing.** Improper handling may cause damage of glass door or injury.
-  **Be careful not to pinch your fingers in the slide rail or the lock mechanism of the drawer.** It may cause injury.

ENVIRONMENTAL CONDITIONS





This equipment is designed to be safe at least under the following conditions (based on the IEC 61010-1):

- Indoor use;
- Altitude up to 2000 m;
- Temperature 5 °C to 40 °C;
- Maximum relative humidity 80 % for temperature up to 31 °C decreasing linearly to 50 % relative humidity at 40 °C;
- Mains supply voltage fluctuations up to ± 10 % of the nominal voltage;
- Transient overvoltages up to the levels of OVERVOLTAGE CATEGORY II;
- Temporary OVERVOLTAGES occurring on the mains supply;
- Applicable pollution degree of the intended environment (POLLUTION DEGREE 2 in most cases);

The above is not a performance specification. Refer to “PERFORMANCE” page for the performance specifications and the environmental conditions.



SYMBOLS ON UNIT

The following symbols are attached to the unit. The table describes the meaning of the symbols.

	This symbol is attached to covers that access high-voltage electrical components to prevent electric shock. Only a qualified engineer or service personnel should be allowed to open these covers.
	This symbol indicates that caution is required. Refer to product documentation for details.
	This symbol indicates incorrect usage could lead to a fire hazard.
	This symbol indicates an earth.
	This symbol means "ON" for a power switch.
○	This symbol means "OFF" for a power switch.

REFRIGERATOR COMPONENTS

Main body

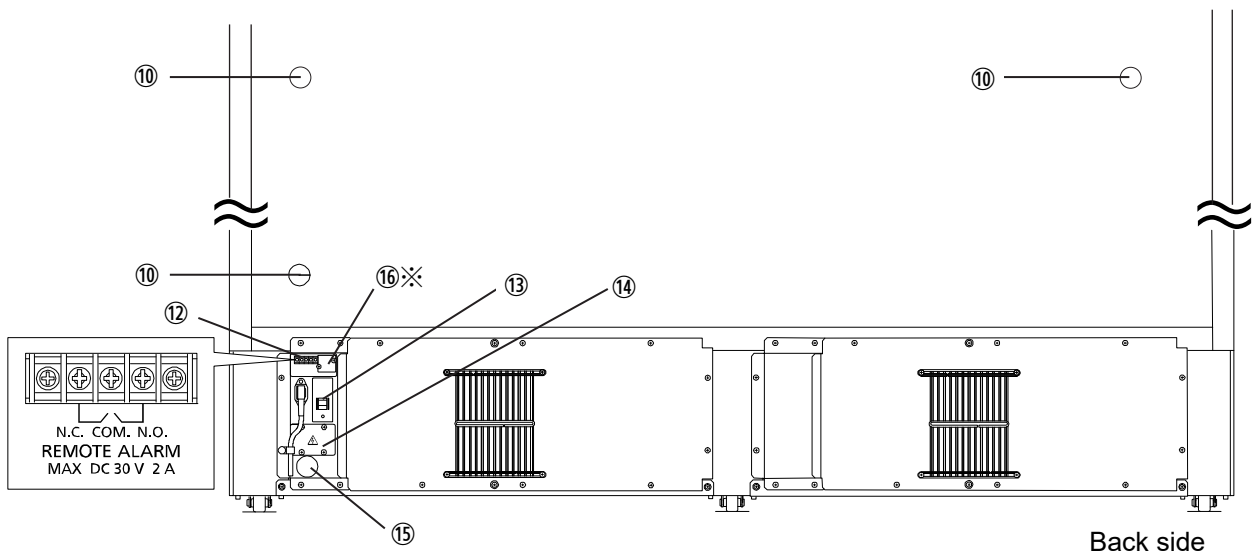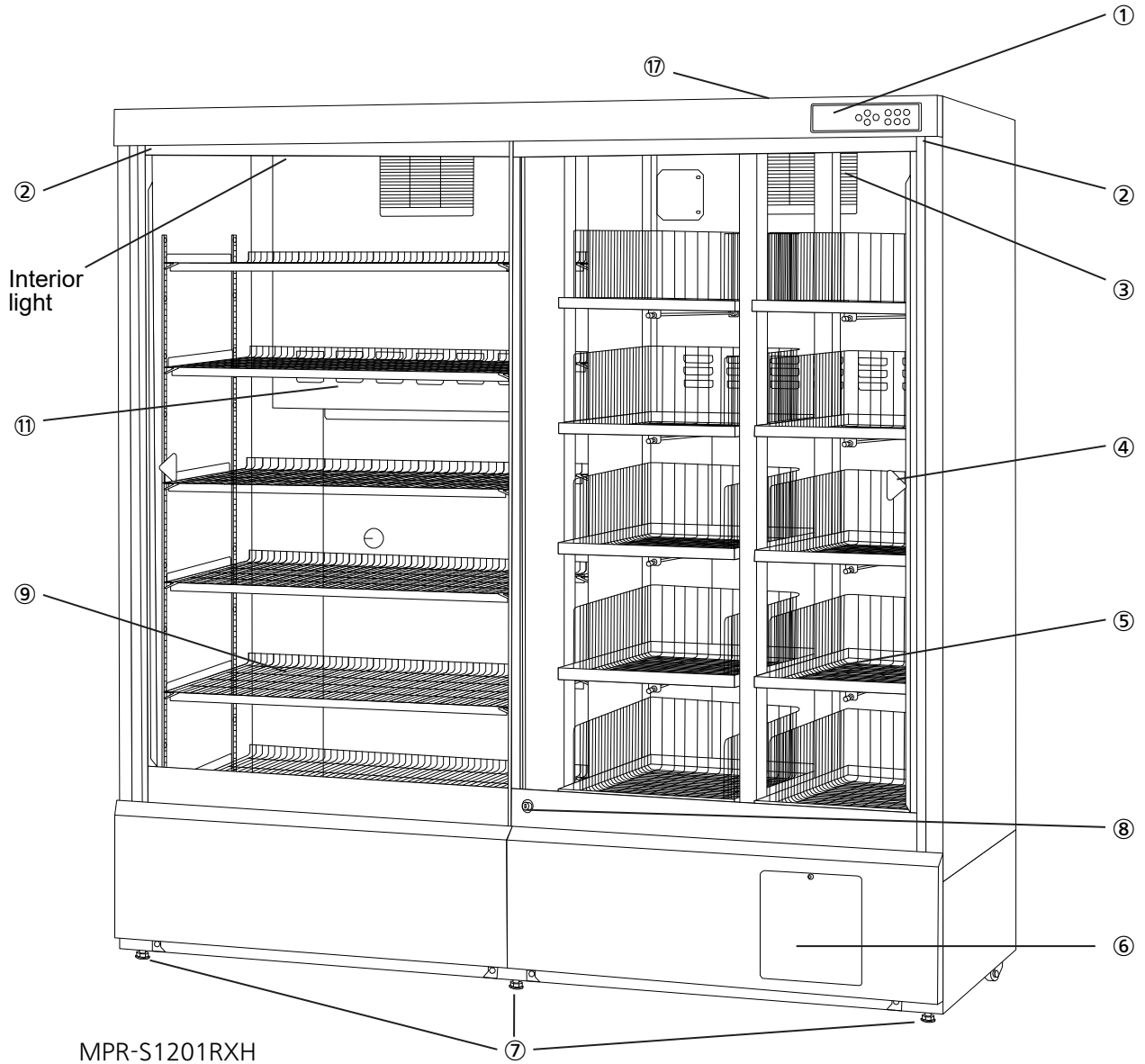


REFRIGERATOR COMPONENTS

- ① **Control panel:** The chamber temperature and other alarms/functions can be set using the keys on the control panel. The operational status can be checked on the temperature display and indicator [page 14].
- ② **Door switch:** This switch detects the door open/close status. The LED lamp on the control panel blinks orange when the door is open.
- ③ **Air intake vent:** This is an air intake vent for circulating the air in the chamber. Do not block this vent.
✦Blocking this vent may cause unstable chamber temperature.
- ④ **Glass door:** Sliding type. The recessed portion on the rail enables the self-closing of the door. Hold the handle indicated by the hand mark to open and close. The glass is double-glazed construction.
✦The condensation may be found on the door frame and on the glass surface depending on the installation environment. This is not a malfunction. This is not unusual.
✦The condensation on the door is dropped on the rail and exhausted to the evaporating tray. However, the water that falls on the rail is not completely drained.
- ⑤ **Drawer (right side of MPR-S1201RXH):** The allowable load is 20 kg per drawer [page 20].
Note: Never pull out all of the multiple drawers at a time. It may cause the unit to tip over.
- ⑥ **Space for temperature recorder:** A temperature recorder (optional) can be mounted here so that the chamber temperature can be recorded automatically [page 61].
✦Contact our sales representative or agent to arrange the installation.
- ⑦ **Levelling foot (front):** These are screw bolts used to install and fix the unit. Adjust the height of the levelling feet by turning the screw bolts until the two front casters are lifted away from the floor [page 17].
- ⑧ **Keyhole:** Push and turn the key clockwise through 90 degree to securely lock the door.
- ⑨ **Shelf:** The location is selectable. Set the shelves horizontally and securely. The allowable load is 50 kg per shelf.
✦Items to be stored in the chamber must be placed on the shelves. Do not put stored items directly on the bottom of the chamber.
- ⑩ **Access port:** This port is used to pass the sensor or cable of measuring equipment, or the sensor of a temperature recorder (optional) to the chamber.
✦Replace the insulation and the rubber caps when the access port is not in use. Improper replacement may cause an increase in chamber temperature or condensation around the access port.
- ⑪ **Air exhaust vent:** This is an air exhaust vent for circulating the air in the chamber. Do not block this vent.
✦Blocking this vent may cause unstable chamber temperature.
- ⑫ **Remote alarm terminal (backside):** A remote alarm device (separately available) can be connected to this terminal. The remote alarm relays the alarm to an operator in a remote location if the unit is unattended [page 53].
✦Contact our sales representative or agent to arrange the installation
- ⑬ **Circuit breaker (backside):** Before starting the refrigerator, check the switch is on. (ON=" | ", OFF="○")
✦The round button under the circuit breaker is a leakage test button. The operation check of the circuit breaker can be performed by pressing this button. But note the power supply to the unit is disconnected when this button is pressed.
- ⑭ **Space for interface board:** An interface board (optional) can be mounted here.
✦Contact our sales representative or agent to arrange the installation.
- ⑮ **Port for servicing (back, left):** This is a access port for servicing. Do not remove the rubber cap.
- ⑯ **Switch for power failure alarm:** This is a switch for power failure alarm [page 61]. Normally, set the switch to ON. Turn off the switch when the refrigerator is not in use.
✦This switch is provided with the optional battery kit for power failure alarm.
- ⑰ **USB port (top):** This port may be used for future enhancements.

REFRIGERATOR COMPONENTS

Control panel

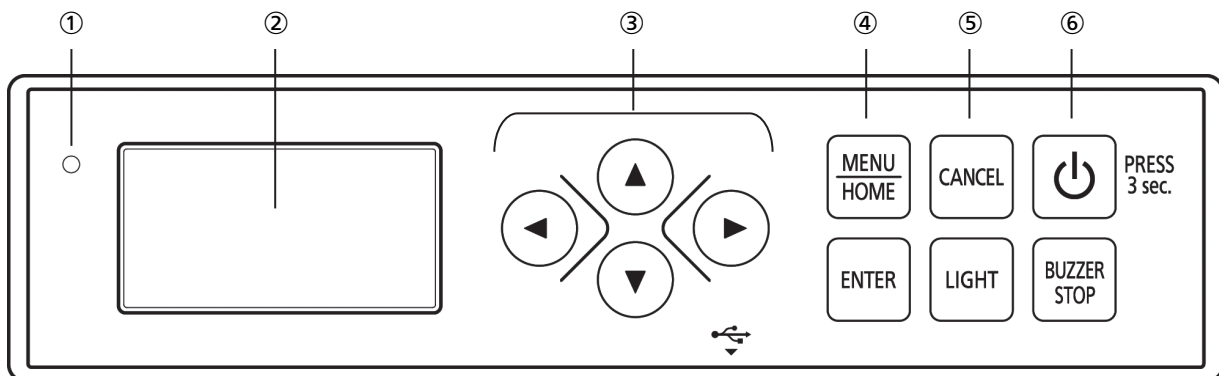

① LED indicator:

In normal operation: The green LED lamp lights up.

During alarm condition: The red LED lamp blinks.

During notification: The orange LED lamp blinks.

During standby status: The orange LED lamp lights up.

② **Temperature display:** The present chamber temperature, log chart, or input screen, etc. are displayed.

③ Up/down/left/right key ($\Delta \nabla \triangleleft \triangleright$):

On the screen "other than the home screen": pressing this key moves the cursor, etc.

④ Menu key (MENU/HOME):

On the "home screen (temperature display screen)": pressing this key opens the menu screen. It is possible to change various setting on the menu screen.

On the "other than the home screen": pressing this key also returns to the home screen.


⑤ **Cancel key (CANCEL):** On the setting screen, pressing this key changes into the screen of one level up.

⑥ **Standby key:** The normal status and standby status is changed by pressing this key for three seconds. The operation stops during standby. Pressing this key for three seconds during normal operation displays the standby screen and makes the LED indicator lamp light up orange. Pressing this key for three seconds during standby status appears the home screen (temperature display) and lights up the green LED lamp.

◇ Under standby status, the refrigerator operation is stopped, and the chamber temperature rises.

◇ The standby key cannot be locked even if the keypad lock is ON [page 26].

REFRIGERATOR COMPONENTS

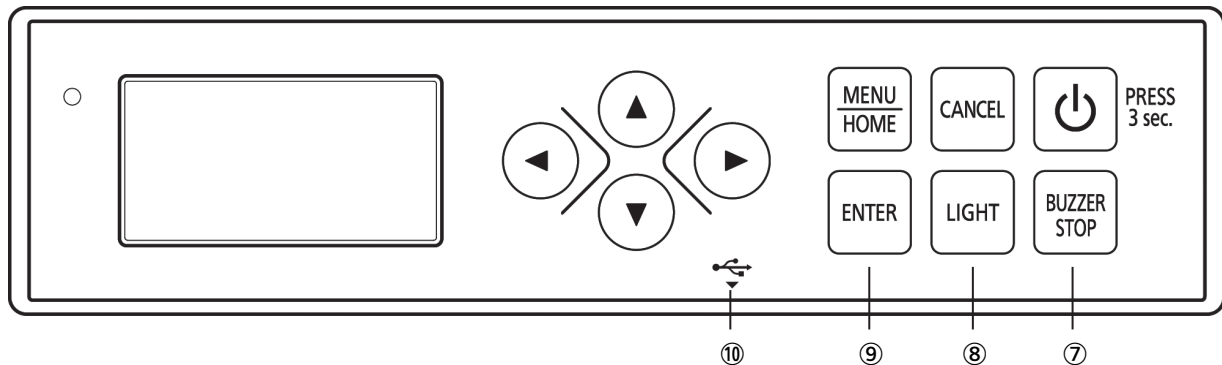

⑦ **Buzzer stop key (BUZZER STOP):** The buzzer is silenced by pressing this key when the buzzer sounds [page 54 and page 55].

✧The alarm status of remote alarm terminal is not cancelled by pressing the buzzer stop key (BUZZER STOP) since the remote alarm is not linked with the buzzer (The buzzer is silenced). Refer to “Setting the remote alarm” on page 45 to link the remote alarm with the buzzer.

✧The buzzer and the remote alarm operate again after certain period (buzzer suspended period) when the same alarm status is continued [page 43].

The chamber temperature is displayed for five seconds by pressing this key during power failure alarm (when battery kit for power failure alarm is installed).

✧Pressing this key for three seconds in home screen activates the alarm test. Pressing this key again finishes the alarm test.

⑧ **Light key (LIGHT):** The light turns on by pressing this key when the light is off. The light turns off by pressing this key when the light is on.

✧This key is invalid when the setting of interior light is ON or OFF.

⑨ **Enter key (ENTER):** Press this to select the menu item. Press this key to enter the required value during the setting procedure.

⑩ **USB port (under side):** Insert a USB flash drive to export data logs.

✧USB flash drives with capacity of 32 GB or less that employ the FAT16/FAT32 file system are supported. USB flash drives that require passwords cannot be used. Do not insert devices other than USB flash drives into the USB port.

INSTALLATION SITE

This unit must be installed in a location which meets all the conditions described below.

◇ If the unit is installed in a location which does not meet the conditions, its specified performance may not be achieved or malfunctions and accidents may occur.

■ A location not exposed to direct sunlight

Avoid any location which is exposed to direct sunlight. Installing the unit in a location exposed to direct sunlight may reduce its cooling performance.

■ A well-ventilated (airy) location

In order to ensure ventilation, leave clearances of at least 10 cm around the unit (at the left, right, top and back). Blocking the ventilation may reduce the unit's cooling performance or cause malfunctions.

■ A location away from sources of heat

Avoid any location which is close to a major source of heat (such as a heater or boiler). Installing the unit near a major source of heat may reduce the unit's cooling performance.

■ A location with minimal changes in temperature

Avoid any location where the ambient temperature is subject to sudden changes. If the unit is installed in a location where the ambient temperature is subject to sudden changes, it will not be possible to achieve a stable cooling performance.

■ A flat surface where the floor is also capable of bearing the total combined weight (product + optional accessories + stored items)

Install the unit on a flat level surface even and which is capable of bearing the total combined weight (product + optional accessories + stored items). If the unit is installed where the surface is uneven or where the unit will be inclined at an angle, the unit will be unstable, and accidents or injuries may occur and/or vibration or unnecessary noise may be generated.

■ A location with minimal humidity

Install the unit in a location where the relative humidity is less than 80 %R.H. Installing the unit in a very humid location may cause earth faults and/or electric shock. If installed at a humidity of 80% or higher, condensation may form on the unit.

■ A location free of flammable or corrosive gases

Avoid any location exposed to flammable or corrosive gases. Flammable or corrosive gases can cause explosions and/or a fire. Furthermore, corrosion of electrical parts may cause insulation to be reduced and result in earth faults and/or electric shock.

■ A location where nothing can fall onto the unit

Avoid locations where objects may fall onto the unit. Objects falling and hitting the unit may cause it to break down or fail.

INSTALLATION

1. Preparations after unpacking

Remove all the tape used to secure the doors and interior parts, and leave the doors open for a short while for ventilation.

If any surfaces of the outer cabinet are dirty, wipe the surface using a cloth moistened with a diluted neutral dish-washing detergent.

✧Using an undiluted solution of detergent may cause the unit's plastic areas to crack. Follow the directions on the detergent for details of dilution.

✧After wiping the unit using the diluted detergent, be absolutely sure to wipe the surfaces with a cloth dipped in clean water to remove traces of the detergent. After this, be absolutely sure to wipe the surfaces with a dry cloth, allowing the surfaces of the outer cabinet to dry out completely, and then proceed with the installation.

2. Securing and levelling the unit using the levelling feet

Rotate the front levelling feet until the casters are raised above the floor surface [Fig. 1].

✧Use a tool (spanner, wrench, etc.) to rotate the levelling foot.

In addition, rotate the levelling feet slightly clockwise or anticlockwise, and adjust them so that the unit is completely level.

✧When the casters are raised from the floor surface, the unit will be secured. If they are left touching the floor, the unit may accidentally move when its door is opened or closed.

3. Securing the unit by using the fixtures



Use the fixtures on the rear panel of the unit, and secure the unit to a wall with a strong rope or chain [Fig. 2].

4. Preventing electric shock by earthing the unit

When installing the unit, be absolutely sure to earth (ground) it. Earthing is necessary to prevent electric shocks resulting from deterioration of electrical insulations.

✧This unit comes with a 3-pin plug having one earth pin. Earthing is not required in the case of a 3-pin power outlet equipped with an earth contact.

✧If the power outlet is not a 3-pin outlet equipped with an earth contact, ask a qualified contractor to do the earthing work.

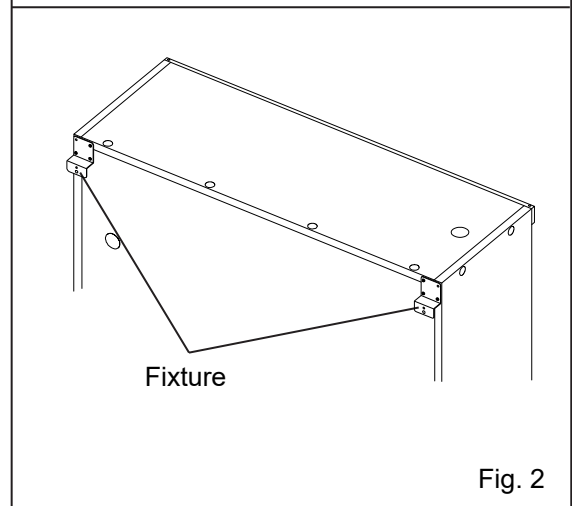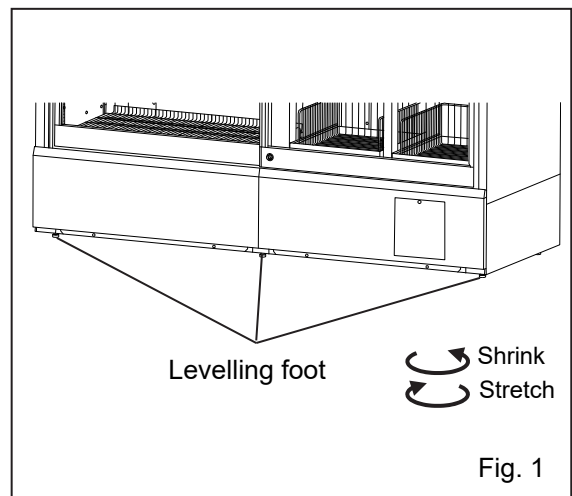

INSTALLATION

5. Setting up the shelves

The shelves are packaged at the bottom of the chamber. Set the shelves firmly in place on the shelf stoppers at the standard locations as marked [Fig. 3].

<Important>

Do not place anything on the bottom of the chamber. Do not put stored items directly on the bottom of the chamber.

How to remove the glass door

Move the glass door to the center position when removing the glass door

✧Take care to handle the glass door carefully after removing. Rough handling may cause damage of the glass door or injury.

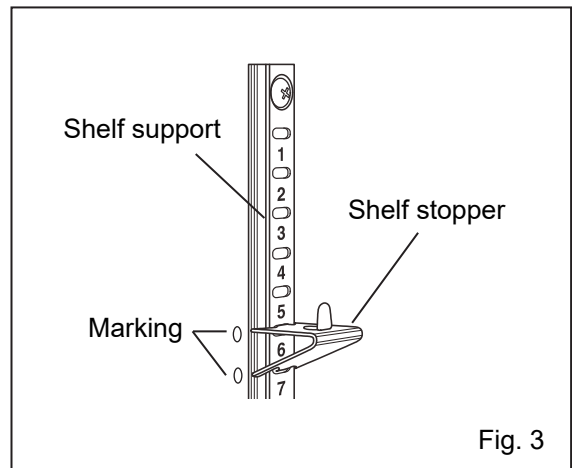


Fig. 3

6. Installing the side guard

To prevent stored items from falling, install the provided side guard on the center side of the shelf by inserting it from above. [Fig. 4].

✧The number of side guard is as follows:

MPR-S1201XH : 12 pcs. (6 already installed on the shelves)

MPR-S1201RXH : 6 pcs. (3 already installed on the shelves)

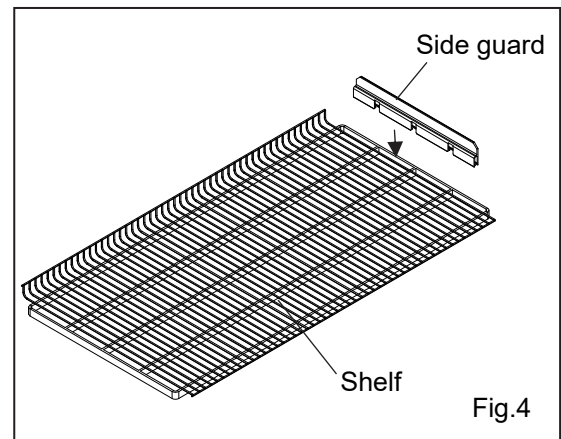


Fig.4

PLACEMENT OF STORED ITEMS

Do not place stored items so as to obstruct the air circulation in the chamber. Obstruction of air circulation can cause the freezing of stored items or can impair the temperature distribution in the chamber.

Air intake vent


Do not block the air intake vent with the stored items. Care should be taken not to put any paper or plastic paper or plastic packaging materials near the air intake vent. These may be sucked in and block the air intake vent.

Air exhaust vent

Do not block the air exhaust by the stored items near the air exhaust vent. Care should be taken not to place stored items near to the air exhaust vent. The stored items may become frozen when the refrigerator is operating at temperature settings lower than 5 °C.

Load line [see figure below]

The storage area is limited to ensure adequate air flow in the chamber. The label "load line" is attached to the right and left sides of the chamber. Always place stored items within the area limited by these labels.

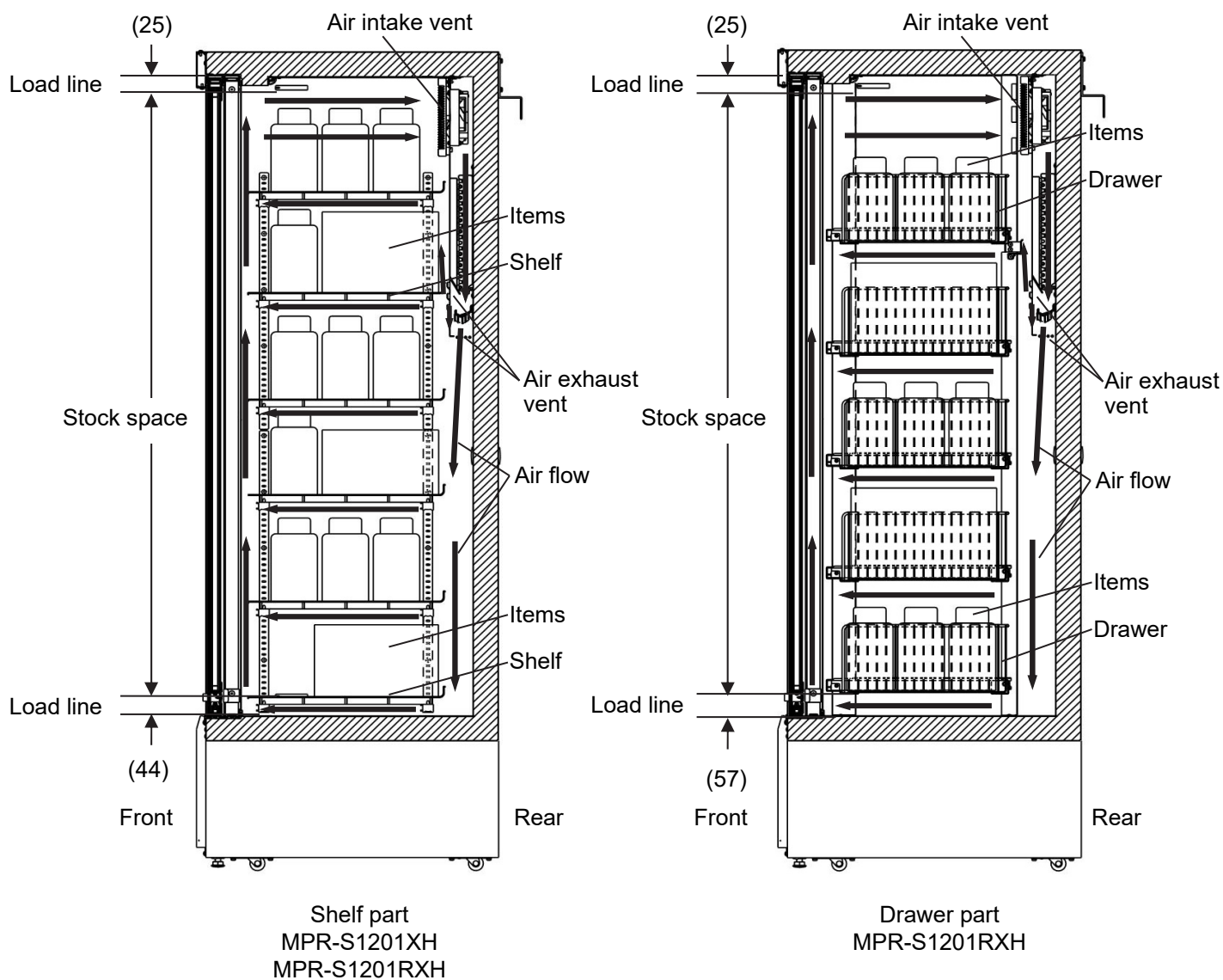

<Side view>

USE OF SHELVES AND DRAWERS

<Shelves>

Do not place anything on the bottom of the chamber. Do not put stored items directly on the bottom of the chamber.

◇Be careful not to drop the stored items from the front of the shelves when placing the stored items on the shelves and when removing the stored items from the shelves.

<Drawers>

Do not place anything on the bottom of the chamber. Do not put stored items directly on the bottom of the chamber.

◇Pay attention to the step on the drawer handle when placing the stored items in the drawer and when removing the stored items from the drawer.

When pulling out a drawer

Pull out a drawer whilst pushing the unlock button [Fig. 1] on the front of the drawer.

When pushing-in a drawer



Push a drawer all the way in until a “click” is heard. The drawer is locked.

When removing a drawer

Pull out a drawer whilst pushing the unlock button [Fig. 1] on the front of the drawer, lift the drawer slightly at the hooked part of the roller [Fig. 2] and take out the drawer.

When installing a drawer

Lift the front of the drawer slightly, insert it into the rail [Fig. 2], and push it in until a “click” is heard. The drawer is locked.

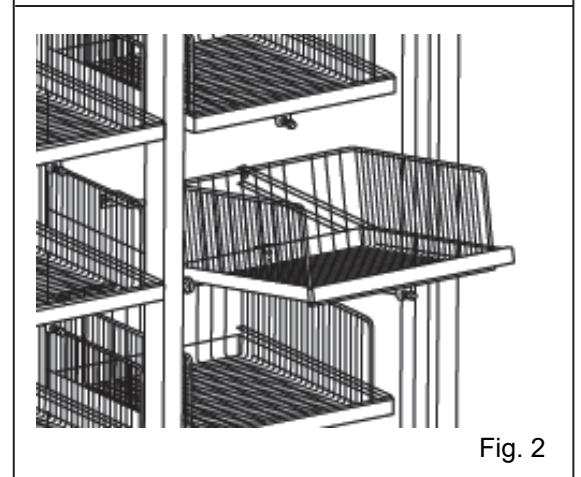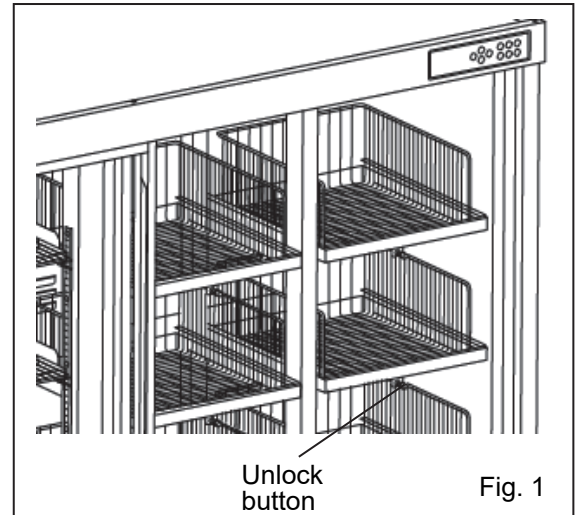

START-UP PROCEDURE

Follow this procedure for the initial operation of the unit and for consequent operations (after temporary stoppage for cleaning, maintenance or moving) .

◇After a power failure, the unit will restart operation automatically with the same settings as before the power failure.

1. Check the circuit breaker switch is ON.

2. Connect the power supply cord to the dedicated power source with the appropriate rating with the chamber empty.

3. Press the standby key on the control panel for more than three seconds.

▶ The home screen (temperature display screen) is displayed. The red LED lamp blinks. (This is caused by the high-temperature alarm function and is not a malfunction.)

4. Turn on the battery switch of the power failure alarm (when the optional battery kit for power failure alarm is installed).

◇“N09:Battery Err” is displayed on the message column when the battery switch of power failure alarm is OFF. The message disappears when the battery switch of power failure alarm is turned ON.

5. Set the date and time displayed on the temperature display to the current date and time [pages 51 to 52].

6. Set the desired chamber temperature [page 25].

◇The factory setting of refrigerator chamber temperature is 5 °C .

7. Open the door to check that the interior light is on.

8. Using the temperature display, check that the chamber temperature has cooled to the set temperature.

9. Gradually place the material inside the chamber.

◇Putting a large amount of material into the chamber at one time causes the temperature to rise.

◇Do not block the air intake vent nor air exhaust vent in the chamber.

10. Set the selectable parameter (keypad lock, alarm temperature, delay of door alarm buzzer, etc.) as required [page 24].

OPERATION DURING POWER FAILURE

This product operates as follows even during a power failure.

	Battery kit for power failure alarm	
	Installed	Not installed
The remote alarm activates.	○	○
The clock continues to operate.	○	○
The power failure alarm (alarm indicator/alarm buzzer) activates.	○*1	×
The chamber temperature is displayed for five seconds by pressing the buzzer stop key.	○*1	×
The log continued to be saved.	○*1	×

*1: When the battery is flat, these functions do not work.

AFTER RECOVERY FROM POWER FAILURE

After recovery from a power failure, the operation will resume automatically with the same settings as before the power failure. There is no need to re-set. However, always check the operational status after recovery.


◇The settings are memorized during power failure by nonvolatile memory.

HOME SCREEN (Temperature display screen)

When the standby key is pressed for more than three seconds, the following home screen (temperature display screen) is displayed.

✦For the details of standby key, refer to item ⑥ on page 14.

✦Following shows the integer display mode.

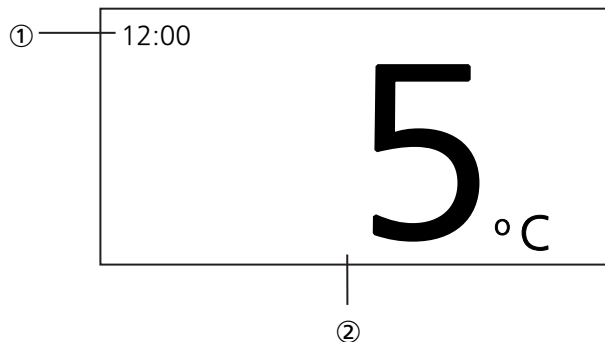

① Message column

Normally, the time (24-hour notation) is displayed. Under alarm status or self-diagnosis function, the error code and message are displayed.

② Temperature display


The current chamber temperature is displayed. During the temperature alarm, the chamber temperature blinks.

<Change of display mode>

On the home screen, the temperature display is selectable; integer display mode or decimal display mode.

Following shows an example of decimal representation mode.

✦For the setting of temperature display, refer to page 39.

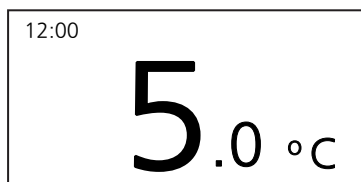

In both the integer display and decimal display, the chamber temperature is displayed repeatedly at certain interval when there is no key operation (sleeping mode).

Pressing any key cancels the repeated display of chamber temperature.

MENU SCREEN

The menu screen is displayed by pressing the menu key (MENU/HOME) while the home screen (temperature display screen) is displayed. Following shows the menu screen.



① Cursor

This is used for selecting the menu. The cursor moves by pressing the up/down key (Δ / ∇).

② Chamber set temperature

Select this menu when changing the refrigerator set temperature. Moving the cursor to this menu and pressing the enter key (ENTER) opens the screen for setting the refrigerator temperature.

③ Keypad lock

Select this menu to set the keypad lock. Moving the cursor to this menu and pressing the enter key (ENTER) opens the screen for setting the keypad lock.

✦The setting can be locked to avoid accidental changes. When the keypad lock is ON, a change of setting is not accepted even if the key on the control panel is operated.

④ Temp. Max / Min

Select this menu to check the maximum or minimum chamber temperature during every 12 hours or 24 hours. Moving the cursor to this menu and pressing the enter key (ENTER) opens the screen for selecting the maximum or minimum temperature display.

⑤ Data Log

Select this menu to display the log on the screen or export the log to a USB flash drive. Moving the cursor to this menu and pressing the enter key (ENTER) opens the “Data Log” screen.

⑥ Alarms & Controls

Select this menu to set the high/low temperature alarm, or other parameters. Moving the cursor to this menu and pressing the enter key (ENTER) opens the “Alarms & controls” screen.

⑦ Temperature display

The setting of chamber temperature is displayed.

MENU SCREEN

Following shows the screen display and function at each screen under Menu level.

Note: The unit will return from setting mode to the home screen automatically after 90 seconds if no key is operated (auto-return function). In this case, the setting is not changed.

Screen	Function	Reference page
"Menu" screen		
"Set Temp." screen		
Set Temp.	Setting the chamber temperature	25
"Keypad Lock" screen		
Keypad Lock ON/OFF	Setting the keypad lock (ON/OFF)	26 to 27
"Temp. Max/Min" screen		
Temp. Max/Min	Display of maximum and minimum chamber temperature	28
Max/Min Reset	Reset of maximum and minimum chamber temperature	29
Max/Min Interval	Setting the interval for maximum/minimum temperature display	30
"Data Log" screen		
Data Log Chart	Chart display of chamber temperature or door operation log	31 to 32
Data Log Export	Exporting the chamber temperature and door operation log	33
Data Log Setting	Setting the log interval, setting of device ID	34 to 35
Event	Display of event log	36
Event Export	Exporting the event log	37
"Alarms & Controls" screen		
1/10 Temp. Display	Setting the decimal display of temperature	39
"Alarm Setting" screen		
High Temp.	Setting the high-temperature alarm	40
Low Temp.	Setting the low-temperature alarm	41
Door Alarm Delay	Setting the delay of door alarm	42
Ring Back Delay	Setting the buzzer suspended period	43
Alarm Volume	Setting the buzzer volume	44
Remote ON/OFF	Setting the remote alarm	45
"Light Setting" screen		
Interior Light	Setting the interior light	46
Lighting Time	Setting the lighting time of the interior light	47
Comp. Delay	Setting the compressor delay	48
Rail Heater	Setting the rail heater operation	49
"Date & Time" screen		
Date Format	Setting the date display (format)	50
Date	Setting the date	51
Time	Setting the time	52

SETTING THE CHAMBER TEMPERATURE

Set the chamber temperature as required to keep the stored material at an appropriate temperature for a long period of time.

- Setting range of chamber temperature: between 2 °C and 14 °C
- Initial setting (factory setting): 5 °C

1. On the home screen, press the menu key (MENU/HOME).

▶ The Menu screen is displayed [Fig. 1].

2. With the cursor on the Set Temp. , press the enter key (ENTER).

▶ The Set Temp. screen is displayed, and the current setting (05) is displayed [Fig. 2].


3. By pressing the right or left key, move the cursor to select the digit to be changed. Then change the figure by pressing the up or down key.

4. Press the enter key (ENTER).


▶ The setting is memorized and the screen is changed to the Menu screen.

5. Press the menu key (MENU/HOME) to display the home screen.

◇ The display returns to the home screen automatically when 90 seconds has passed without any key operation.

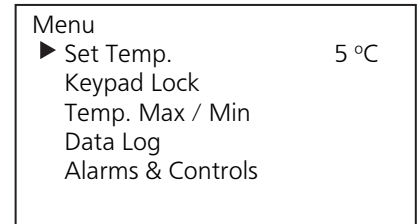

[Fig. 1]

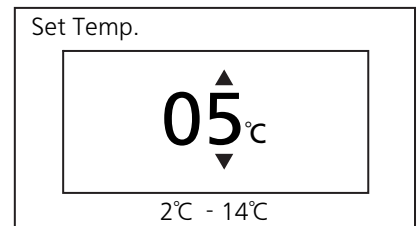

[Fig. 2]


SETTING THE KEYPAD LOCK

The setting can be locked to avoid accidental changes. When the keypad lock is ON, a change of setting is not accepted even if the key on the control panel is operated.


1. On the home screen, press the menu key (MENU/HOME).
 - ▶ The Menu screen is displayed.
2. By pressing the up or down key, move the cursor to the Keypad Lock [Fig. 1] and then press the enter key (ENTER).
 - ▶ The Keypad Lock ON/OFF screen is displayed, and the current setting (OFF) is displayed [Fig. 2].
3. By pressing the up or down key, change the setting to ON.
 - ◇ The setting is changed to ON or OFF when the up or down key is pressed.
4. Press the enter key (ENTER).
 - ▶ The Keypad Lock PW screen is displayed [Fig. 3].
5. Input the password (6 digits) and press the enter key (ENTER).
 - ▶ The Confirm PW screen is displayed [Fig. 4].
6. Again input the password (6 digits) and press the enter key (ENTER).
 - ▶ The display returns to the Menu screen and **LOCK** is displayed on the upper right [Fig. 5].
 - ◇ The number (6 digits) entered here will be the key unlock password.




[Fig. 1]

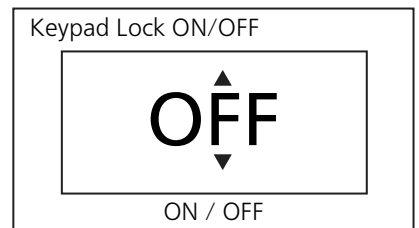

[Fig. 2]

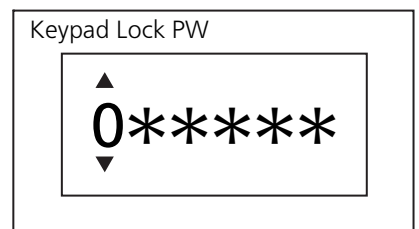


[Fig. 3]

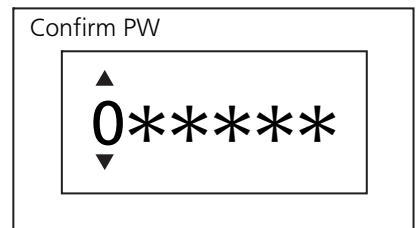


[Fig. 4]

Note: The following screen will be displayed when the second password does not comply with the first password.

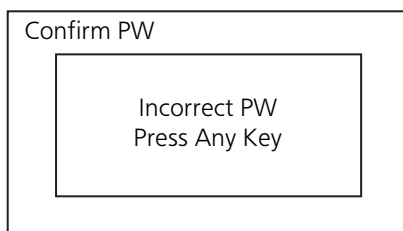

By pressing any key on the control panel, the display returns to the Menu screen. Start from the above procedure 2.

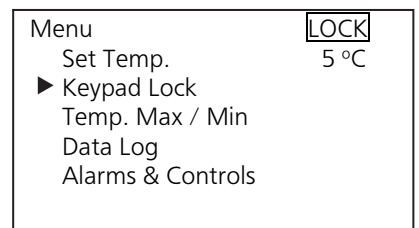

[Fig. 5]


7. Press the menu key (MENU/HOME) to display the home screen.
 - ◇ The display returns to the home screen automatically when 90 seconds has passed without any key operation.

RELEASING THE KEYPAD LOCK

When the keypad lock is OFF, a change of setting is accepted.

1. On the home screen, press the menu key (MENU/HOME).
▶ The Menu screen is displayed.
2. By pressing the up or down key, move the cursor to the Keypad Lock [Fig. 1] and then press the enter key (ENTER).
▶ The Keypad Lock ON/OFF screen is displayed, and the current setting (ON) is displayed [Fig. 2].
3. By pressing the up or down key, change the setting to OFF.
✧ The setting is changed to ON or OFF when the up or down key is pressed.
4. Press the enter key (ENTER).
▶ The Keypad Lock PW screen is displayed [Fig. 3].
5. Input a password (6 digits) and press the enter key (ENTER).
▶ The display returns to the Menu screen and LOCK on the upper right is disappeared [Fig.4].


Note: The following screen will be displayed when the password does not comply with the set password.

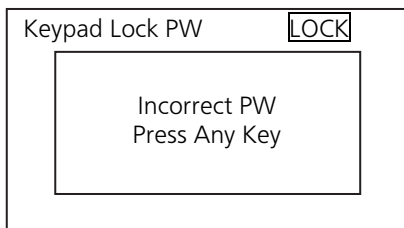


By pressing any key on the control panel, the display returns to the Menu screen. Start from the above procedure 2.

Note: If you forget the password, please contact your sales representative or agent.


6. Press the menu key (MENU/HOME) to display the home screen.
✧ The display returns to the home screen automatically when 90 seconds has passed without any key operation.

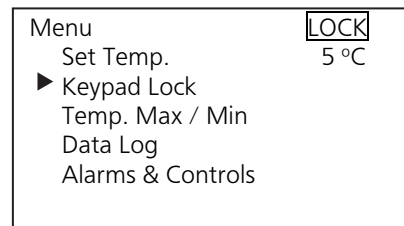

[Fig. 1]

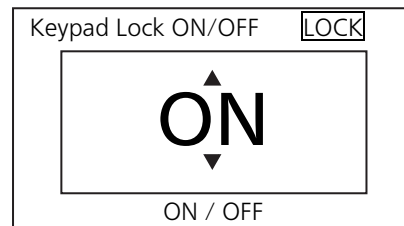

[Fig. 2]

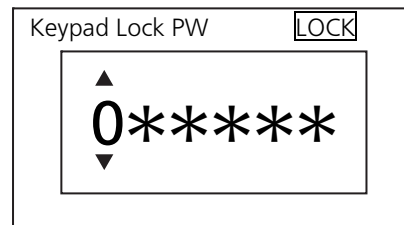

[Fig. 3]



[Fig. 4]


DISPLAY OF MAX/MIN TEMPERATURE

On the temperature display, the maximum and minimum chamber temperature for every 24 or 12 hours can be checked.

- Initial setting (factory setting): 24 hours

1. On the home screen, press the menu key (MENU/HOME).


▶ The Menu screen is displayed.

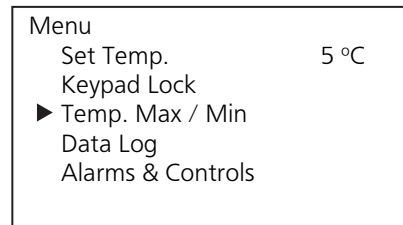

[Fig. 1]

2. By pressing the up or down key, move the cursor to the Temp. Max/Min [Fig. 1] and then press the enter key (ENTER).

▶ The Temp. Max/Min screen is displayed [Fig. 2].

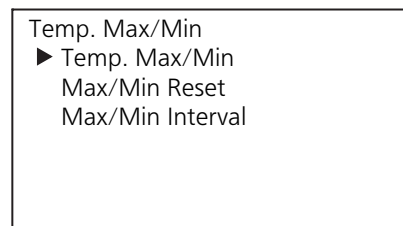


[Fig. 2]

3. With the cursor on the Temp. Max/Min [Fig. 2], press the enter key (ENTER).

▶ The Temp. Max/Min screen is displayed, and the maximum and minimum refrigerator temperature are displayed [Fig. 3].

◇ The display on the top (03/01 12:00) shows the start date and time of the monitor period (for 24 or 12 hours). The start date is March 1 and time is 12:00.

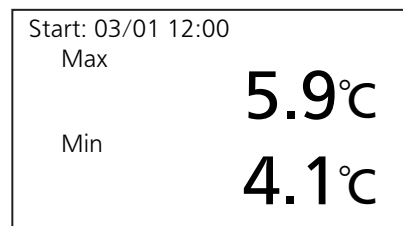

[Fig. 3]


4. Press the menu key (MENU/HOME) to display the home screen.

◇ The display returns to the home screen automatically when 90 seconds has passed without any key operation.


RESET OF MAX/MIN TEMPERATURE

The maximum and minimum chamber temperature can be reset.


1. On the home screen, press the menu key (MENU/HOME).
▶ The Menu screen is displayed.
2. By pressing the up or down key, move the cursor to the Temp. Max/Min [Fig. 1] and then press the enter key (ENTER).
▶ The Temp. Max/Min screen is displayed [Fig. 2].
3. By pressing the up or down key, move the cursor to the Max/Min Reset [Fig. 3] and then press the enter key (ENTER).
▶ The Max/Min Reset screen is displayed, and the current setting (NO) is displayed [Fig. 4].
4. By pressing the up or down key, change the setting to YES.
✧ The setting is changed to NO or YES when the up or down key is pressed.
5. Press the enter key (ENTER).
▶ The setting is memorized and the screen is changed to the Temp. Max/Min screen.
6. Press the menu key (MENU/HOME) to display the home screen.
✧ The display returns to the home screen automatically when 90 seconds has passed without any key operation.

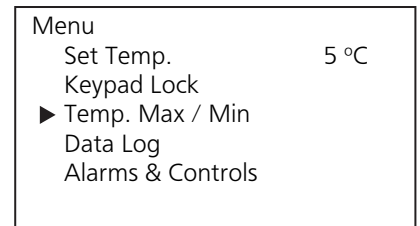


[Fig. 1]

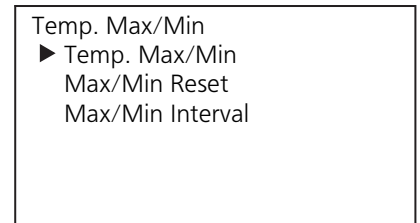

[Fig. 2]

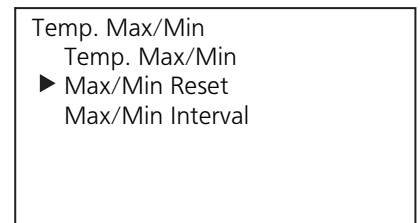

[Fig. 3]

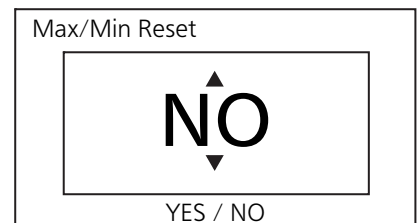

[Fig. 4]

SETTING THE PERIOD FOR MAX/MIN TEMPERATURE

The period for monitoring the maximum and minimum chamber temperature is selectable. The monitor period is 24 hours or 12 hours.


The maximum and minimum chamber temperature for 24 hours are displayed when 24 hours is selected.

The maximum and minimum chamber temperature for 12 hours are displayed when 12 hours is selected.

- Initial setting (factory setting): 24 hours

1. On the home screen, press the menu key (MENU/HOME).


▶ The Menu screen is displayed.

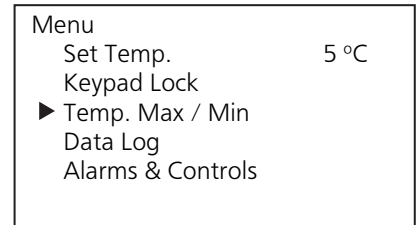

[Fig. 1]

2. By pressing the up or down key, move the cursor to the Temp. Max/Min [Fig. 1] and then press the enter key (ENTER).


▶ The Temp. Max/Min screen is displayed.

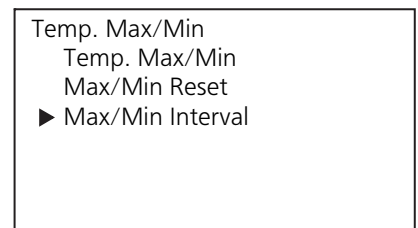

[Fig. 2]

3. By pressing the up or down key, move the cursor to the Max/Min Interval [Fig. 2] and then press the enter key (ENTER).

▶ The Max/Min Interval screen is displayed, and the current setting (24H) is displayed [Fig. 3].

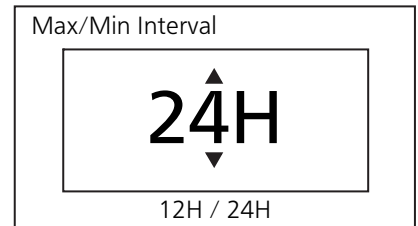

[Fig. 3]

4. By pressing the up or down key, change the setting to 12H.

✧ The setting is changed to 24H or 12H when the up or down key is pressed.

5. Press the enter key (ENTER).

▶ The setting is memorized and the screen is changed to the Temp. Max/Min screen.

6. Press the menu key (MENU/HOME) to display the home screen.

✧ The display returns to the home screen automatically when 90 seconds has passed without any key operation.


DISPLAY AND EXPORT OF DATA LOG

Display & export of temp. and door operation logs

The log of chamber temperature and door operation can be displayed by chart on the screen. Also, those logs can be exported to a USB flash drive. (See page 15 for usable USB flash drives.)

1. On the home screen, press the menu key (MENU/HOME).


▶ The Menu screen is displayed.

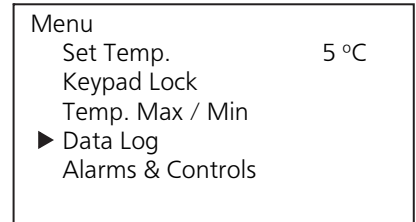

[Fig. 1]

2. By pressing the up or down key, move the cursor to the Data Log [Fig. 1] and then press the enter key (ENTER).


▶ The Data Log screen is displayed [Fig. 2].

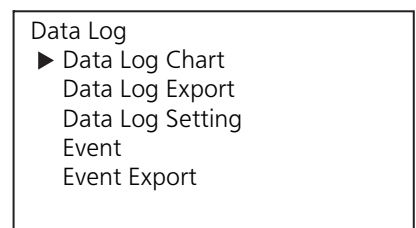

[Fig. 2]

3. With the cursor on the Data Log Chart [Fig. 2], press the enter key (ENTER).

▶ The Temp. screen (chart of refrigerator temperature) is displayed [Fig. 3].

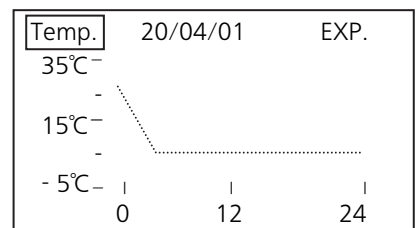


[Fig. 3]

<When displaying the door operation log >

4. With the highlighted Temp. displayed on upper left [Fig. 3], press the enter key (ENTER).

▶ The Data Select screen is displayed [Fig. 4].

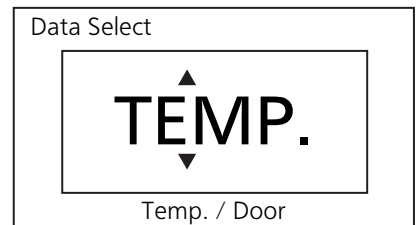

[Fig. 4]

5. By pressing the up or down key, select the data item to be displayed.

◇ The setting is changed to TEMP., or DOOR when the up or down key is pressed.

TEMP.: Log of chamber temperature

DOOR: Log of door operation

6. Press the enter key (ENTER).

▶ The log is displayed in chart.

Note: Fig. 5 shows an example of chart. (OP: door opened, CL: door closed)



[Fig. 5]

<When changing the date for chart display>

7. Press the right key with the data item (Temp. or Door) on upper left is highlighted [Fig. 5].


▶ The date is highlighted.

DISPLAY AND EXPORT OF DATA LOG

8. By pressing the up or down key, select the date of which data is displayed.

◇The date is advanced or moved back by one day when the up or down key is pressed, and the chart of the data for the selected date is displayed.

Note: The Date Select screen is displayed [Fig. 6] when the enter key (ENTER) is pressed instead of the up or down key. Set the date by pressing the up or down key and right or left key.

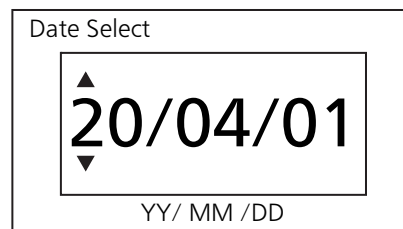

[Fig. 6]

9. Press the enter key (ENTER).

▶ The data for the selected date is displayed in chart.

<When exporting the data for the selected date to a USB flash drive >
(See page 15 for usable USB flash drives.)

10. With the date is highlighted [Fig. 7], press the right key.

▶ EXP. on upper right is highlighted [Fig. 8].

11. Insert a USB flash drive to the USB port.

12. Press the enter key (ENTER).

▶ The logs of chamber temperature and door operation for the selected date are exported with CSV files. During exporting, "Exporting..." is displayed.

Note: Refer to page 59 when a message other than "Exporting..." is displayed.

▶ The chart screen is displayed by pressing the enter key (ENTER).


13. Press any key on the control panel when the screen showing the completion of data exporting is displayed [Fig. 9].

▶ The chart screen is displayed.


14. Remove the USB flash drive from the USB port.

15. Press the menu key (MENU/HOME) to display the home screen.


◇The display returns to the home screen automatically when 90 seconds has passed without any key operation.

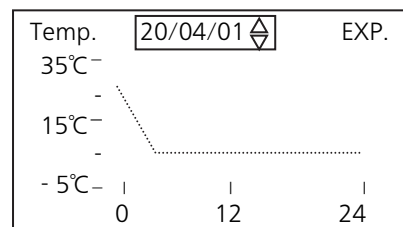

[Fig. 7]

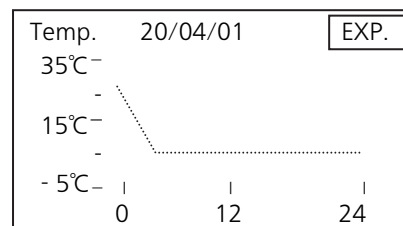

[Fig. 8]

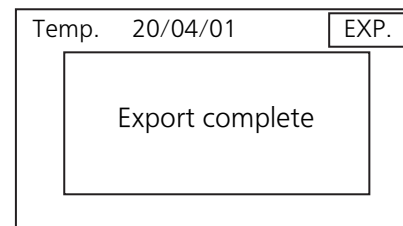

[Fig. 9]

DISPLAY AND EXPORT OF DATA LOG


Export of temperature and door operation logs

All temperature logs and door operation logs saved in a memory or log for a selected date (24 hours) can be exported to a USB flash drive. (See page 15 for usable USB flash drives.)

1. Insert a USB flash drive to the USB port.
2. On the home screen, press the menu key (MENU/HOME).
▶ The Menu screen is displayed.
3. By pressing the up or down key, move the cursor to the Data Log [Fig. 1] and then press the enter key (ENTER).
▶ The Data Log screen is displayed.
4. By pressing the up or down key, move the cursor to the Data Log Export [Fig. 2] and then press the enter key (ENTER).
▶ The Data Log Export screen is displayed [Fig. 3].



[Fig. 1]

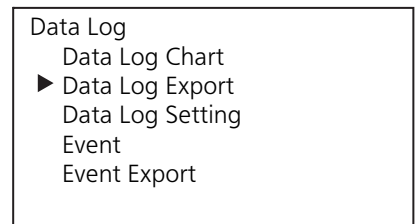

[Fig. 2]

5. With the cursor on the 24 Hours [Fig. 3], press the enter key (ENTER) when the log for a specified date is exported.
▶ The Date Select screen is displayed [Fig.4].




[Fig. 3]

6. By pressing the up, down, right or left key, set a date of which log is exported, and then press the enter key (ENTER).

▶ The log for the set date is exported to the USB flash drive. During exporting, “Exporting...” is displayed.

Note: Refer to page 59 when a message other than “Exporting...” is displayed.

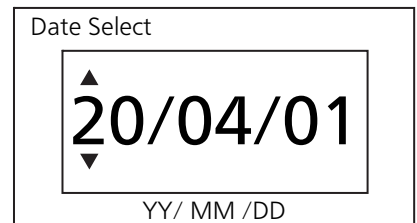

[Fig. 4]

Note: Select All in the above procedure 5 and press the enter key (ENTER) when all memorized logs are exported.

▶ The log folder is created in the USB flash drive, and the exported file is saved in it in CSV format. The exported file name consists of a device ID, date of export and data name. The date of export depends on the setting (refer to page 50).

(e.g.) When exporting data using All (from Jan. 1st, 2019 to Oct. 1st, 2019):

000000_20190101-20191001_DataLog.csv 000000_20190101-20191001_DoorLog.csv

(e.g.) When exporting data using 24 Hours (Jan. 1st, 2019):

000000_20190101_DataLog.csv 000000_20190101_DoorLog.csv

✧The door operation log is exported together.

✧The default setting (000000) is displayed when the device ID is not set.

✧If the file name is duplicated, the data exported earlier is overwritten.

7. “Export complete” is displayed when the export is finished. By pressing the up, down, right or left key, the Data Log Export screen is displayed.

8. Remove the USB flash drive from the USB port.

9. Press the menu key (MENU/HOME) to display the home screen.

✧The display returns to the home screen automatically when 90 seconds has passed without any key operation.

DISPLAY AND EXPORT OF DATA LOG

Setting the log interval

The log interval for the chamber temperature can be set.

◇ When log data reaches the maximum storage amount, the oldest log data is overwritten with the new data.


■ Setting range of log interval: between 1 minute and 15 minutes

■ Initial setting (factory setting): 1 minute (log for approximately 3 months can be stored)

◇ Approximately 3 months of door operation log can be stored. The period may be shortened when the door operation is frequent.

1. On the home screen, press the menu key (MENU/HOME).


▶ The Menu screen is displayed.

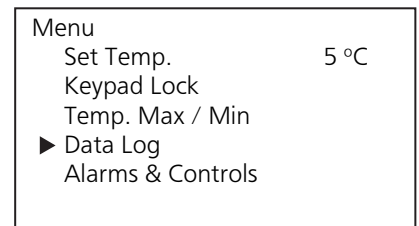

[Fig. 1]

2. By pressing the up or down key, move the cursor to the Data Log [Fig. 1] and then press the enter key (ENTER).


▶ The Data Log screen is displayed.

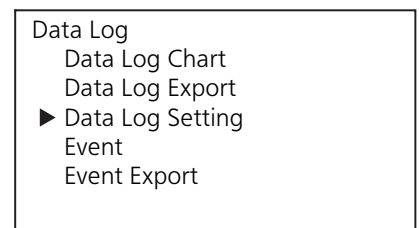

[Fig. 2]

3. By pressing the up or down key, move the cursor to the Data Log Setting [Fig. 2] and then press the enter key (ENTER).


▶ The Data Log Setting screen is displayed [Fig. 3].

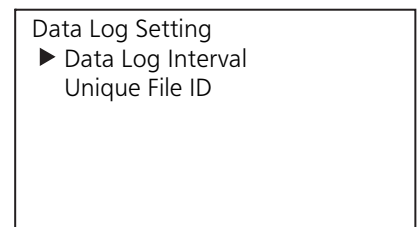

[Fig. 3]

4. With the cursor on the Data Log Interval [Fig. 3], press the enter key (ENTER).

▶ The Data Log Interval screen is displayed [Fig. 4].

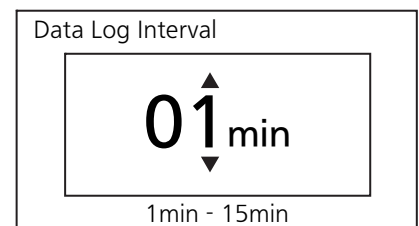

[Fig. 4]

5. By pressing the right or left key, select the digit to be changed and then change the figure by pressing the up or down key.

6. Press the enter key (ENTER).
▶ The setting is memorized and the screen is changed to the Data Log Setting screen.

7. Press the menu key (MENU/HOME) to display the home screen.

◇ The display returns to the home screen automatically when 90 seconds has passed without any key operation.

DISPLAY AND EXPORT OF DATA LOG


Setting the device ID

The unique device ID indicated on the file exported to a USB flash drive (CSV file) can be set.

- Setting range of device ID: 6-digit number and alphabet (capital letter only)
- Initial setting (factory setting): 000000

1. On the home screen, press the menu key (MENU/HOME).

▶ The Menu screen is displayed.

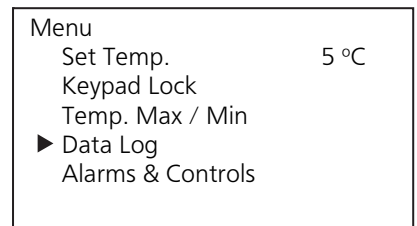

[Fig. 1]

2. By pressing the up or down key, move the cursor to the Data Log [Fig. 1] and then press the enter key (ENTER).

1] and then press the enter key (ENTER).

▶ The Data Log screen is displayed.




[Fig. 2]

3. By pressing the up or down key, move the cursor to the Data Log Setting [Fig. 2] and then press the enter key (ENTER).

▶ The Data Log Setting screen is displayed.

4. By pressing the up or down key, move the cursor to the Unique File ID [Fig. 3] and then press the enter key (ENTER).

▶ The Unique File ID screen is displayed [Fig. 4].

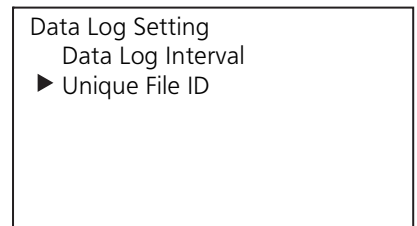


[Fig. 3]

5. By pressing the right or left key, select the digit to be changed and then change the figure or alphabet by pressing the up or down key.

6. Press the enter key (ENTER).

▶ The setting is memorized and the screen is changed to the Data Log Setting screen.

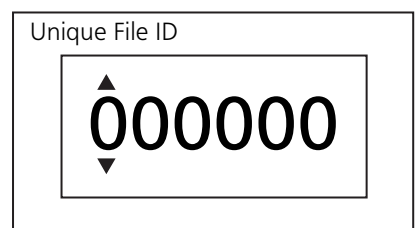

[Fig. 4]


7. Press the menu key (MENU/HOME) to display the home screen.

◇ The display returns to the home screen automatically when 90 seconds has passed without any key operation.

DISPLAY OF EVENT LOG

The event log (max. about 40,000 records) can be displayed on the screen. Following shows the format of the indication.

◇ When the number of logs exceeds the maximum number, the oldest log data is overwritten with the new data.

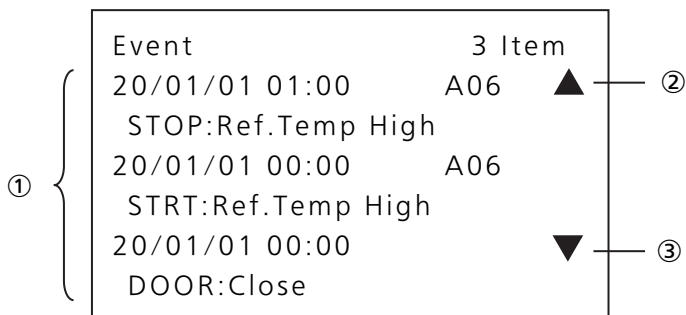

① Display area for event log

One event log is displayed in two lines and three event logs can be displayed on the screen. The upper line shows; a date and time when an event occurred and an error code; a date and time when an event finished and an error code. The lower line shows an error message.

◇ For details of error codes and error messages, refer to page 54 and page 55.

② Up scroll mark (When the number of events is 4 or more)

◇ By pressing the up key, the later event log is displayed when the up scroll mark is displayed.

③ Down scroll mark (When the number of events is 4 or more)

◇ By pressing the down key, the previous event log is displayed when the down scroll mark is displayed.

1. On the home screen, press the menu key (MENU/HOME).


▶ The Menu screen is displayed.



[Fig. 1]

2. By pressing the up or down key, move the cursor to the Data Log [Fig. 1] and then press the enter key (ENTER).


▶ The Data Log screen is displayed [Fig. 2].

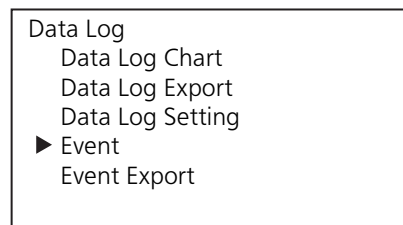

[Fig. 2]

3. By pressing the up or down key, move the cursor to the Event [Fig. 2] and then press the enter key (ENTER).

▶ The Event screen is displayed [Fig. 3].

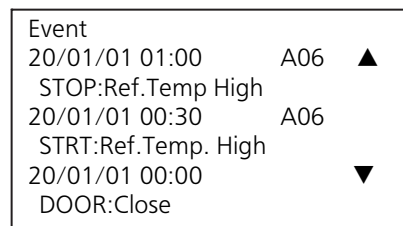

[Fig.3]


4. Press the menu key (MENU/HOME) to display the home screen.

◇ The display returns to the home screen automatically when 90 seconds has passed without any key operation.


EXPORT OF EVENT LOG

The event log for the selected date or all event logs saved in the memory can be exported to a USB flash drive in CSV file. (See page 15 for usable USB flash drives.)

1. Insert a USB flash drive to the USB port.
2. On the home screen, press the menu key (MENU/HOME).
▶ The Menu screen is displayed.
3. By pressing the up or down key, move the cursor to the Data Log [Fig. 1] and then press the enter key (ENTER).
▶ The Data Log screen is displayed.
4. By pressing the up or down key, move the cursor to the Event Export [Fig. 2] and then press the enter key (ENTER).
▶ The Event Export screen is displayed [Fig. 3].
5. With the cursor on the 24 Hours [Fig. 3], press the enter key (ENTER) when the event log for a specified date is exported.
▶ The Date Select screen is displayed [Fig.4].

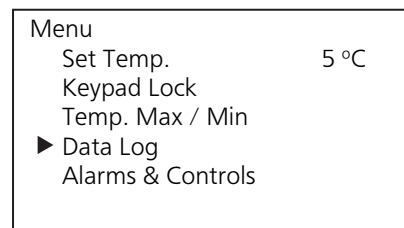

[Fig. 1]

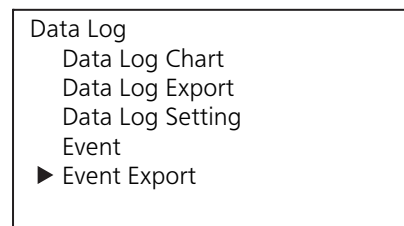


[Fig. 2]



[Fig. 3]

6. By pressing the up, down, right or left key, set a date of which event log is exported, and then press the enter key (ENTER).
▶ The event log for the set date is exported to the USB flash drive. During exporting, “Exporting....” is displayed.
Note: Refer to page 59 when a message other than “Exporting...” is displayed.

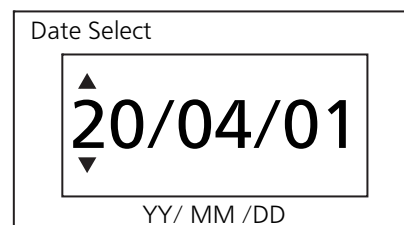

[Fig. 4]

Note: Select All in the above procedure 5 and press the enter key (ENTER) when all memorized event logs are exported.

7. “Export complete” is displayed when the export is finished. By pressing the up, down, right or left key, the Event Export screen is displayed.

▶ The log folder is created in the USB flash drive, and the exported file is saved in it in CSV format. The exported file name consists of a device ID, date of export and data name. The date of export depends on the setting (refer to page 50).

(e.g.) When exporting data using All (from Jan. 1st, 2019 to Oct. 1st, 2019):
000000_20190101-20191001_EventLog.csv

(e.g.) When exporting data using 24 Hours (Jan. 1st, 2019):
000000_20190101_EventLog.csv

◇ The default setting (000000) is displayed when the device ID is not set.

◇ If the file name is duplicated, the data exported earlier is overwritten.

8. Remove the USB flash drive from the USB port.


9. Press the menu key (MENU/HOME) to display the home screen.

◇ The display returns to the home screen automatically when 90 seconds has passed without any key operation.

SETTING THE SELECTABLE PARAMETERS

Selectable parameters

The setting of some parameters (including the alarm setting) can be changed. Following shows the selectable parameters. Refer to the following pages for the details.

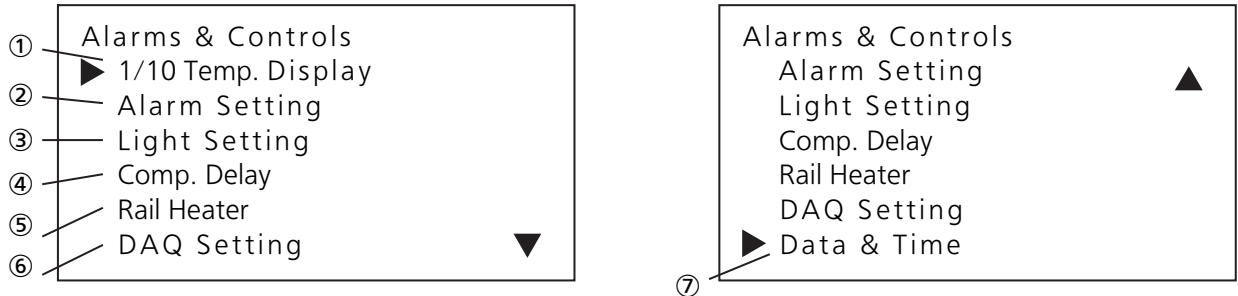

① 1/10 Temp. Display

The temperature display format on the home screen is selectable; integer display or decimal display.

② Alarm Setting

The selectable parameters; high temperature alarm and low temperature alarm, delay of door alarm buzzer, buzzer suspended period, buzzer volume, remote alarm.

③ Interior Light

The setting of interior light; linked with the door operation, always ON or always OFF and setting of lighting period.

④ Comp. Delay

The delay time of compressor start after recovery from the power failure is selectable.

⑤ Rail Heater

The setting of rail heater; Auto change, always OFF or always ON.

⑥ DAQ Setting

The selectable parameters; device ID for data acquisition, communication speed of data acquisition, remote control.

⑦ Date & Time

The selectable parameters; setting of date and time, and format of date/time indication “YY/MM/DD” or “DD/MM/YY”


SETTING THE SELECTABLE PARAMETERS

Setting the temperature display


1. On the home screen, press the menu key (MENU/HOME).
▶ The Menu screen is displayed.
2. By pressing the up or down key, move the cursor to the Alarms & Controls [Fig. 1] and then press the enter key (ENTER).
▶ The Alarms & Controls screen is displayed [Fig. 2].
3. With the cursor on the 1/10 Temp. Display [Fig. 2], press the enter key (ENTER).
▶ The 1/10 Temp. Display screen is displayed [Fig.3].
4. The setting is changed to the integer display (00°C) or decimal display (00.0°C) when the up or down key is pressed.
5. Press the enter key (ENTER).
▶ The setting is memorized and the screen is changed to the Alarms & Controls screen.
6. Press the menu key (MENU/HOME) to display the home screen.
◇ The display returns to the home screen automatically when 90 seconds has passed without any key operation.



[Fig. 1]

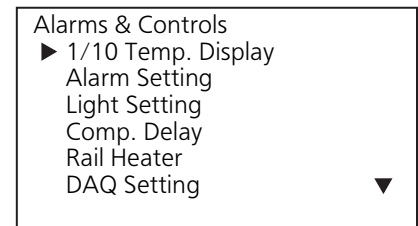

[Fig. 2]

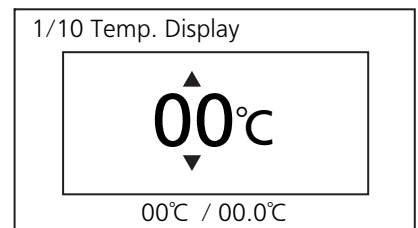

[Fig. 3]

SETTING THE SELECTABLE PARAMETERS

Setting the high-temperature alarm

An abnormality (chamber temperature rise) is indicated by the blinking of the red LED lamp and the chamber temperature display and the buzzer sounding (15 minutes after blinking) if the chamber temperature exceeds the set value of the high-temperature alarm. Always set an appropriate temperature according to the contents to protect stored items from damage resulting from an increase in temperature.

- Setting range for high-temperature alarm:
between 2 °C and 14 °C higher than the chamber set temperature
- Initial setting (factory setting): 5 °C higher than the chamber set temperature

1. On the home screen, press the menu key (MENU/HOME).


▶ The Menu screen is displayed.



[Fig. 1]

2. By pressing the up or down key, move the cursor to the Alarms & Controls [Fig. 1] and then press the enter key (ENTER).


▶ The Alarms & Controls screen is displayed.

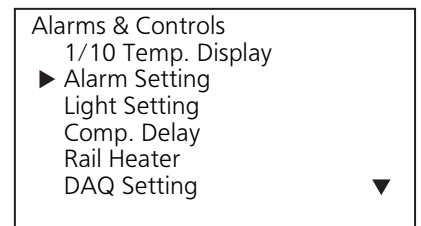

[Fig. 2]

3. By pressing the up or down key, move the cursor to the Alarm Setting [Fig. 2] and then press the enter key (ENTER).


▶ The Alarm Setting screen is displayed.

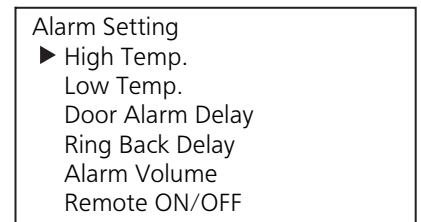

[Fig. 3]

4. With the cursor on the High Temp. [Fig. 3], press the enter key (ENTER).

▶ The High Temp. screen is displayed, and the current setting (05) is displayed [Fig. 4].

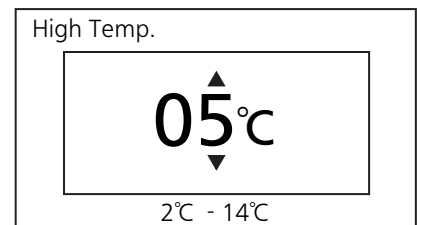

[Fig. 4]

5. By pressing the right or left key, select the digit to be changed and then change the figure by pressing the up or down key.

6. Press the enter key (ENTER).

▶ The setting is memorized and the screen is changed to the Alarm Setting screen.

7. Press the menu key (MENU/HOME) to display the home screen.

◇ The display returns to the home screen automatically when 90 seconds has passed without any key operation.

SETTING THE SELECTABLE PARAMETERS


Setting the low-temperature alarm

An abnormality (chamber temperature drop) is indicated by the blinking of red LED lamp and the chamber temperature display and the buzzer sounding (15 minutes after blinking) if the chamber temperature falls below the set value of low-temperature alarm. Always set an appropriate temperature according to the contents to protect stored items from damage resulting from a drop in temperature.

- Setting range for low-temperature alarm:
between 14 °C and 2 °C lower than the chamber set temperature
- Initial setting (factory setting): 5 °C lower than the chamber set temperature

1. On the home screen, press the menu key (MENU/HOME).


▶ The Menu screen is displayed.

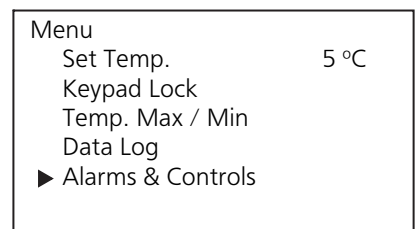

[Fig. 1]

2. By pressing the up or down key, move the cursor to the Alarms & Controls [Fig. 1] and then press the enter key (ENTER).


▶ The Alarms & Controls screen is displayed.

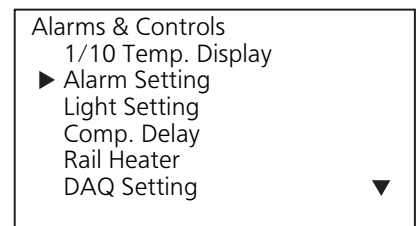

[Fig. 2]

3. By pressing the up or down key, move the cursor to the Alarm Setting [Fig. 2] and then press the enter key (ENTER).


▶ The Alarm Setting screen is displayed.

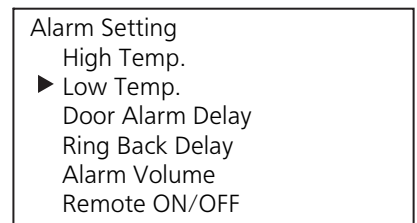

[Fig. 3]

4. By pressing the up or down key, move the cursor to the Low Temp. [Fig. 3] and then press the enter key (ENTER).

▶ The Low Temp. screen is displayed, and the current setting (-05) is displayed [Fig. 4].

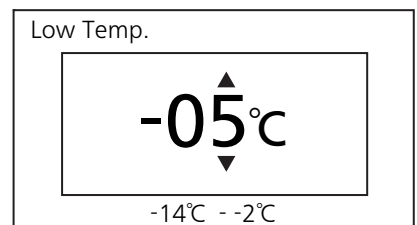

[Fig. 4]

5. By pressing the right or left key, select the digit to be changed and then change the figure by pressing the up or down key.

6. Press the enter key (ENTER).

▶ The setting is memorized and the screen is changed to the Alarm Setting screen.

7. Press the menu key (MENU/HOME) to display the home screen.

◇ The display returns to the home screen automatically when 90 seconds has passed without any key operation.

SETTING THE SELECTABLE PARAMETERS


Setting the door buzzer delay

The red LED lamp blinks and “N02:Door Delay ON” is displayed when the door is opened. After a delay, the display is changed to “N01:Door Open” and the buzzer sounds to indicate that the door is open. The delay time (between the blinking of the red LED lamp and the activation of the buzzer) can be changed. Set an appropriate delay time according to the condition of use to prevent the rise of chamber temperature occurring as a result of the door being left open for too long.

- Setting range of delay time: between 0 and 15 minutes (0 means no delay)
- Initial setting (factory setting): 2 minutes

1. On the home screen, press the menu key (MENU/HOME).


▶ The Menu screen is displayed.

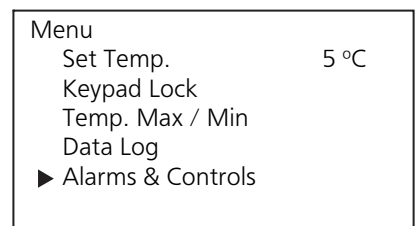

[Fig. 1]

2. By pressing the up or down key, move the cursor to the Alarms & Controls [Fig. 1] and then press the enter key (ENTER).


▶ The Alarms & Controls screen is displayed.

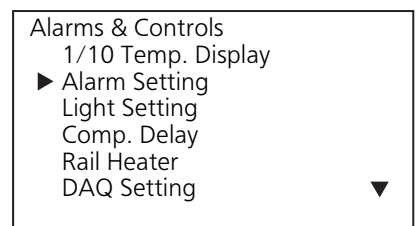

[Fig. 2]

3. By pressing the up or down key, move the cursor to the Alarm Setting [Fig. 2] and then press the enter key (ENTER).


▶ The Alarm Setting screen is displayed.

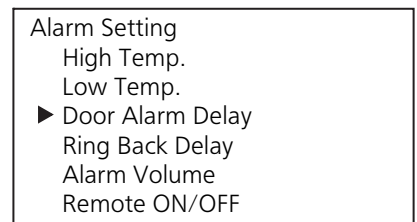

[Fig. 3]

4. By pressing the up or down key, move the cursor to the Door Alarm Delay [Fig. 3], press the enter key (ENTER).

▶ The Door Alarm Delay screen is displayed, and the current setting (02) is displayed [Fig. 4].

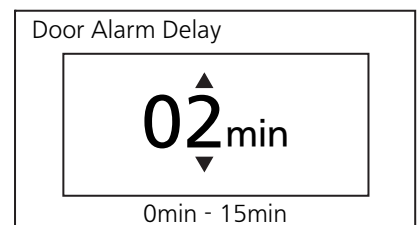

[Fig. 4]

5. By pressing the right or left key, select the digit to be changed and then change the figure by pressing the up or down key.

6. Press the enter key (ENTER).

▶ The setting is memorized and the screen is changed to the Alarm Setting screen.

7. Press the menu key (MENU/HOME) to display the home screen.

◇ The display returns to the home screen automatically when 90 seconds has passed without any key operation.

SETTING THE SELECTABLE PARAMETERS

Setting the buzzer suspended period

An abnormality is indicated again by the buzzer sounding after a certain period has elapsed (the buzzer recovery time) if the alarm mode continues even after the buzzer has already been silenced by pressing the buzzer stop key (BUZZER STOP). Always set the “buzzer suspended period” to avoid incorrect recognition of alarm mode.


- Setting range of buzzer suspended period: between 0 and 60 minutes in 10-minute intervals
- Initial setting (factory setting): 30 minutes

◇The buzzer does not sound again when the setting is 0 (but if the other alarm is occurred, the buzzer sounds). The setting should be between 10 and 60 to ensure the safety of stored items.

◇The setting during the alarm status is invalid even if the buzzer is stopped. It is valid when the alarm status is canceled or the buzzer is sounds resulting from a new alarm mentioned on page 54 and page 55.

1. On the home screen, press the menu key (MENU/HOME).


▶ The Menu screen is displayed.

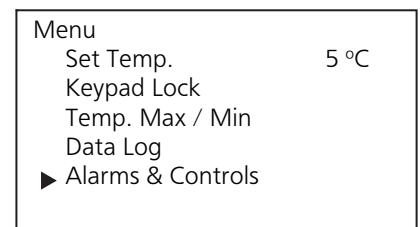

[Fig. 1]

2. By pressing the up or down key, move the cursor to the Alarms & Controls [Fig. 1] and then press the enter key (ENTER).


▶ The Alarms & Controls screen is displayed.

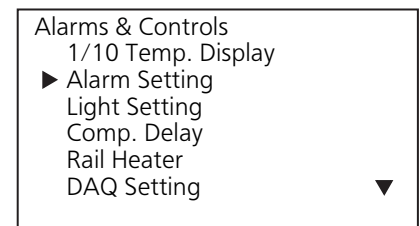

[Fig. 2]

3. By pressing the up or down key, move the cursor to the Alarm Setting [Fig. 2] and then press the enter key (ENTER).


▶ The Alarm Setting screen is displayed.

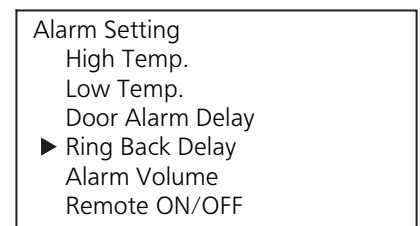

[Fig. 3]

4. By pressing the up or down key, move the cursor to the Ring Back Delay [Fig. 3] and then press the enter key (ENTER).

▶ The Ring Back Delay screen is displayed, and the current setting (30) is displayed [Fig. 4].

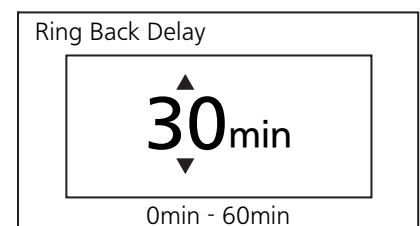

[Fig. 4]

5. By pressing the up or down key, change the figure.

6. Press the enter key (ENTER).

▶ The setting is memorized and the screen is changed to the Alarm Setting screen.

7. Press the menu key (MENU/HOME) to display the home screen.

◇The display returns to the home screen automatically when 90 seconds has passed without any key operation.

SETTING THE SELECTABLE PARAMETERS


Setting the buzzer volume

The buzzer volume is selectable. Select the appropriate volume taking into consideration of the environment of installation site.

- Setting range of buzzer volume: Low (small = standard), High (loud)
- Initial setting (factory setting): Low (small = standard)

1. On the home screen, press the menu key (MENU/HOME).


▶ The Menu screen is displayed.

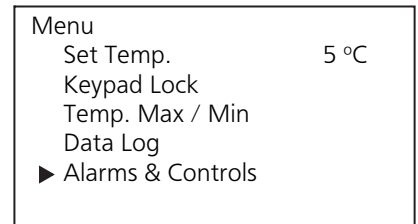

[Fig. 1]

2. By pressing the up or down key, move the cursor to the Alarms & Controls [Fig. 1] and then press the enter key (ENTER).


▶ The Alarms & Controls screen is displayed.

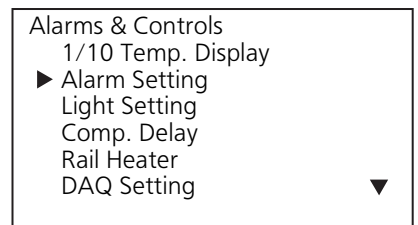

[Fig. 2]

3. By pressing the up or down key, move the cursor to the Alarm Setting [Fig. 2] and then press the enter key (ENTER).


▶ The Alarm Setting screen is displayed.

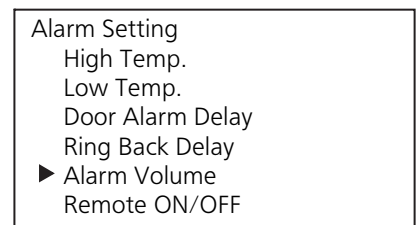

[Fig. 3]

4. By pressing the up or down key, move the cursor to the Alarm Volume [Fig. 3] and then press the enter key (ENTER).

▶ The Alarm Volume screen is displayed, and the current setting (LOW) is displayed [Fig. 4].

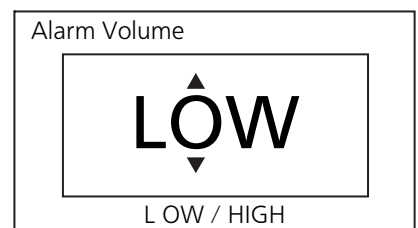

[Fig. 4]

5. By pressing the up or down key, change the setting to HIGH.

◇ The setting is changed to LOW or HIGH when the up or down key is pressed.

6. Press the enter key (ENTER).

▶ The setting is memorized and the screen is changed to the Alarm Setting screen.

7. Press the menu key (MENU/HOME) to display the home screen.

◇ The display returns to the home screen automatically when 90 seconds has passed without any key operation.

SETTING THE SELECTABLE PARAMETERS

Setting the remote alarm

The alarm of this unit can be relayed to a remote location from this unit by connecting the external alarm device (commercial goods) to the remote alarm terminals. It is selectable whether the operation of the external alarm device is linked with the buzzer or not.

In the case of link: The alarm of the external device is cancelled when the buzzer stop key (BUZZER STOP) is pressed during buzzer operation. But the alarm is not cancelled in the case of power failure alarm.

In the case of no-link: The alarm of the external device is not cancelled when the buzzer stop key (BUZZER STOP) is pressed during buzzer operation.

- Setting range of remote alarm: OFF (no-link), ON (link)
- Initial setting (factory setting): OFF (no-link)

1. On the home screen, press the menu key (MENU/HOME).


▶ The Menu screen is displayed.



[Fig. 1]

2. By pressing the up or down key, move the cursor to the Alarms & Controls [Fig. 1] and then press the enter key (ENTER).


▶ The Alarms & Controls screen is displayed.

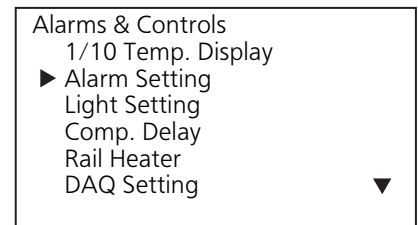

[Fig. 2]

3. By pressing the up or down key, move the cursor to the Alarm Setting [Fig. 2] and then press the enter key (ENTER).


▶ The Alarm Setting screen is displayed.

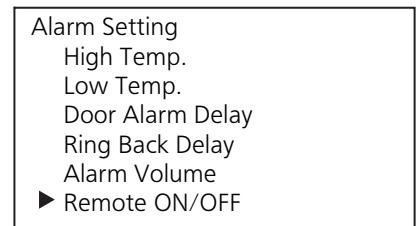

[Fig. 3]

4. By pressing the up or down key, move the cursor to the Remote ON/OFF [Fig. 3] and then press the enter key (ENTER).

▶ The Remote ON/OFF screen is displayed, and the current setting (OFF) is displayed [Fig. 4].

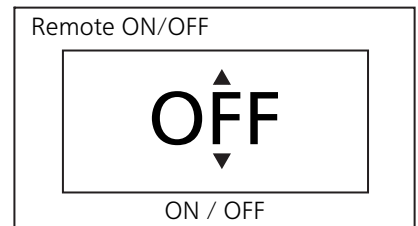

[Fig. 4]

5. By pressing the up or down key, change the setting to ON.

◇ The setting is changed to ON or OFF when the up or down key is pressed.

6. Press the enter key (ENTER).

▶ The setting is memorized and the screen is changed to the Alarm Setting screen.

7. Press the menu key (MENU/HOME) to display the home screen.

◇ The display returns to the home screen automatically when 90 seconds has passed without any key operation.

SETTING THE SELECTABLE PARAMETERS


Setting the interior light

The condition of the interior light can be selected from three options; link with door operation, always ON, always OFF.

- Setting range of interior light: link with door operation (AUTO), always ON (ON), always OFF (OFF)
- Initial setting (factory setting): link with door operation (AUTO)

1. On the home screen, press the menu key (MENU/HOME).


▶ The Menu screen is displayed.

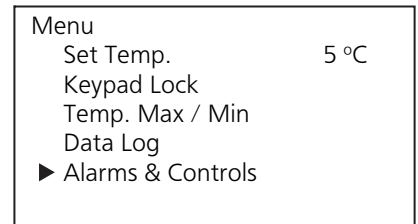

[Fig. 1]

2. By pressing the up or down key, move the cursor to the Alarms & Controls [Fig. 1] and then press the enter key (ENTER).


▶ The Alarms & Controls screen is displayed.

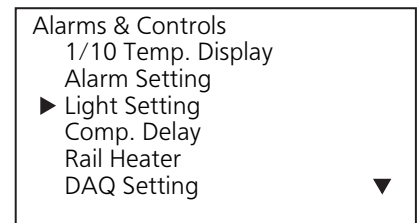

[Fig. 2]

3. By pressing the up or down key, move the cursor to the Light Setting [Fig. 2] and then press the enter key (ENTER).


▶ The Light Setting screen is displayed.

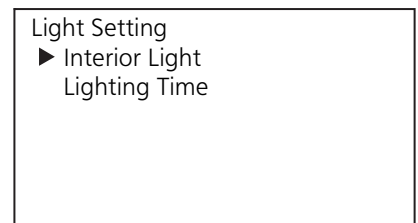

[Fig. 3]

4. With the cursor on the Interior Light [Fig. 3], press the enter key (ENTER).

▶ The Interior Light screen is displayed, and the current setting (AUTO) is displayed [Fig. 4].

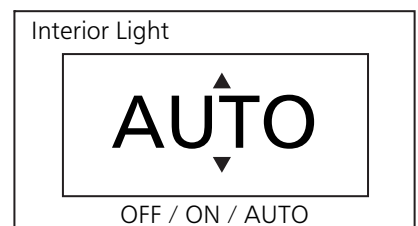

[Fig. 4]

5. By pressing the up or down key, change the setting.

◇ The setting is changed to OFF, ON or AUTO when the up or down key is pressed.

6. Press the enter key (ENTER).

▶ The setting is memorized and the screen is changed to the Alarms & Controls screen.

7. Press the menu key (MENU/HOME) to display the home screen.

◇ The display returns to the home screen automatically when 90 seconds has passed without any key operation.

SETTING THE SELECTABLE PARAMETERS

Setting the lighting time


The lighting time of the interior light can be selected. The interior light turns off automatically when the setting time has passed.

◇The setting is valid only when the setting of interior light is "AUTO".

- Setting range of lighting time: between 1 minute and 15 minutes
- Initial setting (factory setting): 10 minutes

1. On the home screen, press the menu key (MENU/HOME).


▶ The Menu screen is displayed.

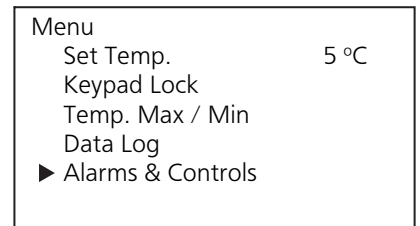

[Fig. 1]

2. By pressing the up or down key, move the cursor to the Alarms & Controls [Fig. 1] and then press the enter key (ENTER).


▶ The Alarms & Controls screen is displayed.

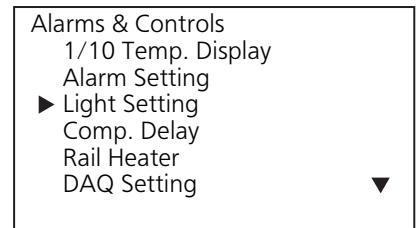

[Fig. 2]

3. By pressing the up or down key, move the cursor to the Light Setting [Fig. 2] and then press the enter key (ENTER).


▶ The Light Setting screen is displayed.

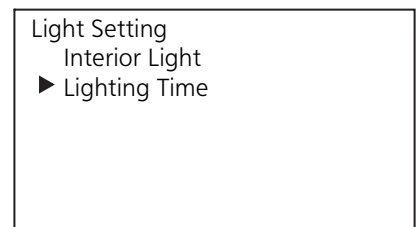

[Fig. 3]

4. By pressing the up or down key, move the cursor to the Lighting Time [Fig. 3] and then press the enter key (ENTER).

▶ The Lighting Time screen is displayed and the current setting (10) is displayed [Fig. 4].

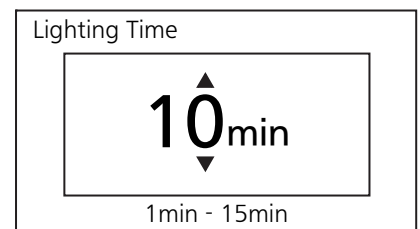

[Fig. 4]

5. By pressing the right or left key, select the digit to be changed and then change the figure by pressing the up or down key.

6. Press the enter key (ENTER).

▶ The setting is memorized and the screen is changed to the Light Setting screen.

7. Press the menu key (MENU/HOME) to display the home screen.

◇The display returns to the home screen automatically when 90 seconds has passed without any key operation.

SETTING THE SELECTABLE PARAMETERS

Setting the compressor delay

The delay time before the compressor starts can be changed to reduce the load on the power line and to facilitate the start-up (reset) of the refrigerator during recovery after power failure.


<Important>

Change the delay time when the multiple devices are connected to one power source.

- Setting range of delay time: between 5 minutes and 15 minutes
- Initial setting (factory setting): 5 minutes

1. On the home screen, press the menu key (MENU/HOME).


▶ The Menu screen is displayed.

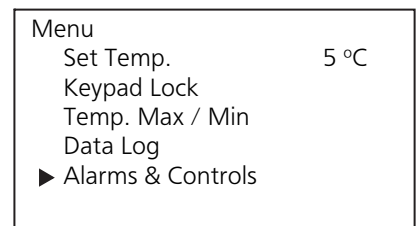

[Fig. 1]

2. By pressing the up or down key, move the cursor to the Alarms & Controls [Fig. 1] and then press the enter key (ENTER).


▶ The Alarms & Controls screen is displayed.

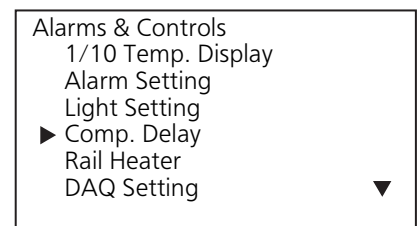

[Fig. 2]

3. By pressing the up or down key, move the cursor to the Comp. Delay [Fig. 2] and then press the enter key (ENTER).

▶ The Comp. Delay screen is displayed, and the current setting (05) is displayed [Fig. 3].

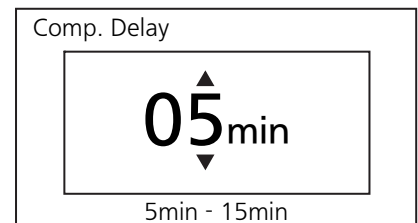

[Fig. 3]

4. By pressing the right or left key, select the digit to be changed and then change the figure by pressing the up or down key.

5. Press the enter key (ENTER).

▶ The setting is memorized and the screen is changed to the Alarms & Controls screen.

6. Press the menu key (MENU/HOME) to display the home screen.

◇ The display returns to the home screen automatically when 90 seconds has passed without any key operation.

SETTING THE SELECTABLE PARAMETERS

Setting the rail heater

This unit has a rail heater installed inside the rail above the door to control condensation on the rail section.

The operation mode can be selected from three options; Auto change, always OFF, and always ON.

◇Auto change automatically controls the heater according to the chamber temperature.

◇Power consumption can be reduced by turning it off all the time, for example, when the product is in an environment where condensation does not occur so often.

- Setting range of operation mode: Auto change (AUTO), always ON (ON), always OFF (OFF)
- Initial setting (factory setting): Auto change (AUTO)

1. On the home screen, press the menu key (MENU/HOME).


▶ The Menu screen is displayed.



[Fig. 1]

2. By pressing the up or down key, move the cursor to the Alarms & Controls [Fig. 1] and then press the enter key (ENTER).

▶ The Alarms & Controls screen is displayed.

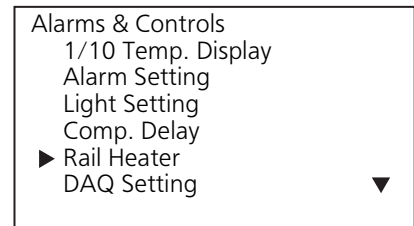


[Fig. 2]

3. By pressing the up or down key, move the cursor to the Rail Heater [Fig. 2] and then press the enter key (ENTER).

▶ The Rail Heater screen is displayed. [Fig. 3].

4. By pressing the up or down key, change the setting.

◇The setting is changed to OFF, ON or AUTO when the up or down key is pressed.

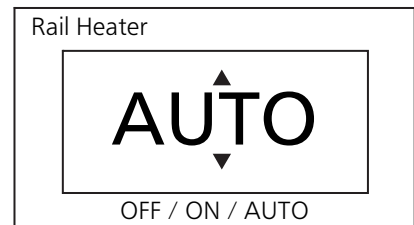

[Fig. 3]

5. Press the enter key (ENTER).

▶ The setting is memorized and the screen is changed to the Alarms & Controls screen.

6. Press the menu key (MENU/HOME) to display the home screen.

◇The display returns to the home screen automatically when 90 seconds has passed without any key operation.

.

SETTING THE SELECTABLE PARAMETERS


Setting the date display (format)

The format of date display is selectable between two types.

- Setting range: year / month / day (YY/MM/DD), or day / month / year (DD/MM/YY)
- Initial setting (factory setting): year / month / day (YY/MM/DD)

1. On the home screen, press the menu key (MENU/HOME).


▶ The Menu screen is displayed.

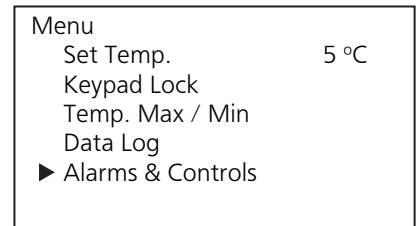

[Fig. 1]

2. By pressing the up or down key, move the cursor to the Alarms & Controls [Fig. 1] and then press the enter key (ENTER).


▶ The Alarms & Controls screen is displayed.

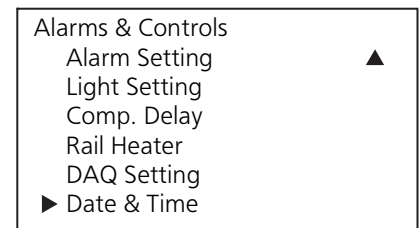

[Fig. 2]

3. By pressing the up or down key, move the cursor to the Date & Time [Fig. 2] and then press the enter key (ENTER).


▶ The Date & Time screen is displayed [Fig. 3].

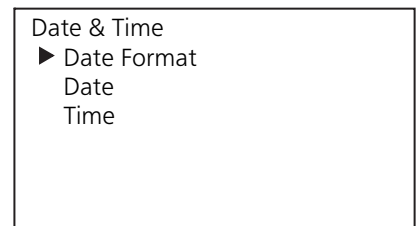

[Fig. 3]

4. With the cursor on the Date Format [Fig. 3], press the enter key (ENTER).

▶ The Date Format screen is displayed, and the current setting (YY/MM/DD) is displayed [Fig. 4].

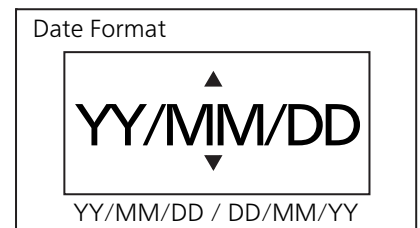

[Fig. 4]

5. By pressing the up or down key, change the setting to DD/MM/YY.

◇ The setting is changed to YY/MM/DD, DD/MM/YY when the up or down key is pressed.

6. Press the enter key (ENTER).

▶ The setting is memorized and the screen is changed to the Date & Time screen.


7. Press the menu key (MENU/HOME) to display the home screen.

◇ The display returns to the home screen automatically when 90 seconds has passed without any key operation.


SETTING THE SELECTABLE PARAMETERS

Setting the date


1. On the home screen, press the menu key (MENU/HOME).
▶ The Menu screen is displayed.
2. By pressing the up or down key, move the cursor to the Alarms & Controls [Fig. 1] and then press the enter key (ENTER).
▶ The Alarms & Controls screen is displayed.
3. By pressing the up or down key, move the cursor to the Date & Time [Fig. 2] and then press the enter key (ENTER).
▶ The Date & Time screen is displayed.
4. By pressing the up or down key, move the cursor to the Date [Fig. 3] and then press the enter key (ENTER).
▶ The Date screen is displayed, and the current date (YY/MM/DD) is displayed [Fig. 4].
5. By pressing the right or left key, select the digit to be changed and then change the figure by pressing the up or down key.
The figure can be changed same as above when the date format is DD/MM/YY.
6. Press the enter key (ENTER).
▶ The setting is memorized and the screen is changed to the Date & Time screen.
7. Press the menu key (MENU/HOME) to display the home screen.
◇ The display returns to the home screen automatically when 90 seconds has passed without any key operation.

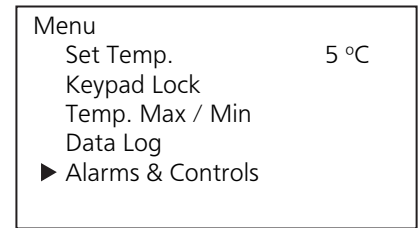


[Fig. 1]

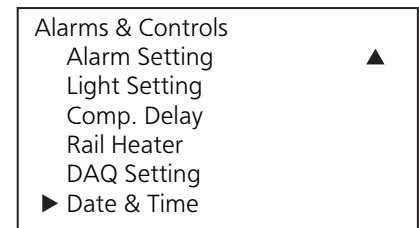

[Fig. 2]

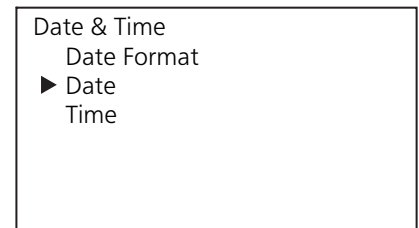

[Fig. 3]

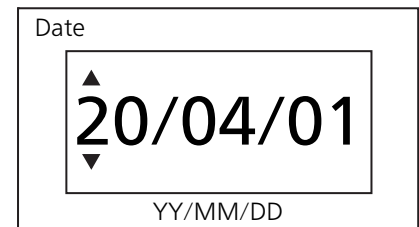

[Fig. 4]


SETTING THE SELECTABLE PARAMETERS

Setting the time

It is recommended to adjust the time periodically since the deviation of about one minutes per a month may be occurred.

1. On the home screen, press the menu key (MENU/HOME).


▶ The Menu screen is displayed.

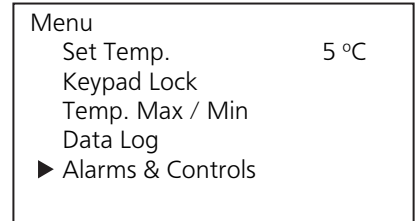

[Fig. 1]

2. By pressing the up or down key, move the cursor to the Alarms & Controls [Fig. 1] and then press the enter key (ENTER).


▶ The Alarms & Controls screen is displayed.

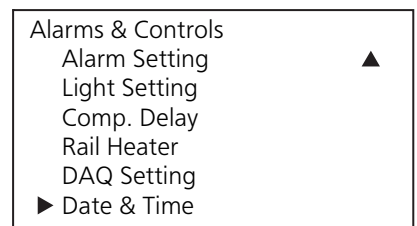

[Fig. 2]

3. By pressing the up or down key, move the cursor to the Date & Time [Fig. 2] and then press the enter key (ENTER).


▶ The Date & Time screen is displayed.

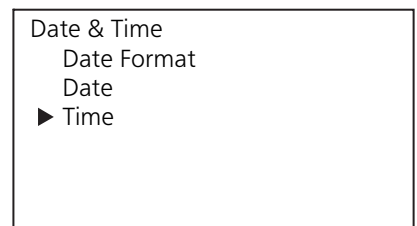

[Fig. 3]

4. By pressing the up or down key, move the cursor to the Time [Fig. 3] and then press the enter key (ENTER).

▶ The Time screen is displayed, and the current time is displayed [Fig. 4].

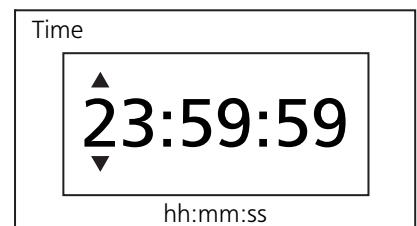

[Fig. 4]

5. By pressing the right or left key, select the digit to be changed and then change the figure by pressing the up or down key.

◇hh:mm:ss means the hour, minute and second can be set in 2-digit.

hh = hour (between 00 and 23), mm = minute (between 00 and 59), ss = second (between 00 and 59)

6. Press the enter key (ENTER).

▶ The setting is memorized and the screen is changed to the Date & Time screen.

7. Press the menu key (MENU/HOME) to display the home screen.

◇The display returns to the home screen automatically when 90 seconds has passed without any key operation.

REMOTE ALARM TERMINAL

The alarm is relayed to a remote location when a remote alarm device (commercially available) is connected to the remote alarm terminal. Installation of a remote alarm device is recommended when the unit is installed in an unattended location so that the operator of the unit is notified of the alarm.

◇Contact our representative or agent to arrange the installation of a remote alarm device.

- Location of remote alarm terminal: left bottom (back side)

- Allowable contact capacity: DC 30 V, 2 A

◇It is recommended to use interface cables with a maximum length of 30 meters.

Contact output

Terminal	At normal	At abnormal
Between COM. and N.O.	open	close
Between COM. and N.C.	close	open

<Important>

The remote alarm terminal will be in alarm mode when the power supply cord is unplugged or the circuit breaker is turned off because this will be interpreted as a power failure.

ALARMS AND SELF-DIAGNOSIS

This unit has the alarm functions and self-diagnosis functions listed below.

An error code and error message are displayed on the message column when an alarm or self-diagnosis occurs.

Contact our sales representative or agent if the alarm mode continues, since the failure of the unit is supposed.

◇Before calling for service, take some precautions to protect the stored items (for example by transferring the items to another refrigerator).

The first letter of the error code means:

A: Alarm - The cooling performance has been significantly degraded. Thus, the chamber temperature may get considerably higher. *1

N: Notification - Cooling performance may have been degraded or there is a possibility of failure other than the cooling performance.

Error code	Error message	Alarm content	LED indicator	Buzzer	Chamber temp. display	Remote alarm *2
A01	Ref.SNSR R Op.	Disconnection of thermal sensor R	Blinking of red lamp	Intermittent tone	ON	Alarm mode
A02	Ref.SNSR R Sh.	Short-circuit of thermal Sensor R	Blinking of red lamp	Intermittent tone	ON	Alarm mode
A03	Ref.SNSR L Op.	Disconnection of thermal sensor L	Blinking of red lamp	Intermittent tone	ON	Alarm mode
A04	Ref.SNSR L Sh.	Short-circuit of thermal Sensor L	Blinking of red lamp	Intermittent tone	ON	Alarm mode
A05	Freeze Warning	Alarm of chamber temp.; equal or lower than 0 °C	Blinking of red lamp	Intermittent tone	Blinking	Alarm mode
A06	Ref.Temp.High	High-temperature alarm	Blinking of red lamp	Intermittent tone	Blinking	Alarm mode
A07	Ref.Temp.Low	Low-temperature alarm	Blinking of red lamp	Intermittent tone	Blinking	Alarm mode
A08*3	Power Failure	Power failure alarm*4	Blinking of red lamp	Intermittent tone	OFF	Alarm mode
A09	INT Fan R Lock	Chamber fan R lock	Blinking of red lamp	Intermittent tone	ON	Alarm mode
A10	INT Fan L Lock	Chamber fan L lock	Blinking of red lamp	Intermittent tone	ON	Alarm mode
A11	CMP Fan R Lock	Compressor fan R lock	Blinking of red lamp	Intermittent tone	ON	Alarm mode
A12	CMP Fan L Lock	Compressor fan L lock	Blinking of red lamp	Intermittent tone	ON	Alarm mode

ALARMS AND SELF-DIAGNOSIS

Error code	Error message	Alarm content	LED indicator	Buzzer	Chamber temp. display	Remote alarm ^{*2}
N01	Door Open	Door alarm	Blinking of orange lamp	Intermittent tone	ON	---
N02	Door Delay ON	Delayed status of door alarm	Blinking of orange lamp	---	ON	---
N03	Ref.Temp.High	High-temperature alarm	Blinking of orange lamp	---	Blinking	---
N04	Ref.Temp.Low	Low-temperature alarm	Blinking of orange lamp	---	Blinking	---
N05	Def.SNSR R Op.	Disconnection of defrost sensor R	Blinking of orange lamp	Intermittent tone	ON	---
N06	Def.SNSR R Sh.	Short-circuit of defrost sensor R	Blinking of orange lamp	Intermittent tone	ON	---
N07	Def.SNSR L Op.	Disconnection of defrost sensor L	Blinking of orange lamp	Intermittent tone	ON	---
N08	Def.SNSR L Sh.	Short-circuit of defrost sensor L	Blinking of orange lamp	Intermittent tone	ON	---
N09 ^{*3}	Battery Err.	Battery error	Blinking of orange lamp	---	ON	---
N10 ^{*3}	Replace Battery	The time for battery replacement	Blinking of orange lamp	---	ON	---
N11	COMM Err.	Communication error	Blinking of orange lamp	---	ON	---
N12	Cool.R Degrade	Degrade of cooling circuit R	Blinking of orange lamp	---	---	---
N13	Cool.L Degrade	Degrade of cooling circuit L	Blinking of orange lamp	---	---	---

*1: The buzzer operate again after certain period when the same alarm status is continued [page 43].

*2: The alarm status is not cancelled by pressing the buzzer stop key (BUZZER STOP) in initial setting (factory setting). Refer to "Setting the remote alarm" on page 45 to link the remote alarm with the buzzer.

*3: When the optional battery kit for power failure alarm is installed.N09 is displayed by turning off the battery switch. N09 may be displayed when the battery is discharged.

*4: The chamber temperature is displayed for five seconds and the buzzer stops when the buzzer stop key (BUZZER STOP) is pressed during the power failure alarm. The red LED lamp continues to blink.

ROUTINE MAINTENANCE

Cleaning of exterior, interior, door rail (including drain port) and accessories

Use a dry cloth to wipe off small amounts of dirt on the outside, inside, door rail (including drain port) of the unit and all accessories. If the outside panels are dirty, clean them with a diluted neutral dishwashing detergent.

Wipe off the condensation on the glass or exterior of the cabinet with a dry soft cloth.

◇Undiluted detergent can damage the plastic components. For the dilution, refer to the instruction of the detergent.

◇After the cleaning with the diluted detergent, always wipe it off with a wet cloth. Then wipe off the cabinet or accessories with a dry cloth.

<Important>

Never pour water onto or into the unit. This may cause electric shock or failure.


Do not use a brush, an acid, a thinner, laundry soap, a powder detergent, or boiling water for cleaning. These may cause damage to painted surfaces or cause perishing of plastic and rubber components. Moreover, do not wipe plastic and rubber components with a volatile material.

Drain port

The drain port is for draining the condensed water that has accumulated on the door rail. The condensed water may not drain when the drain port clogs, and the floor will get wet. Therefore, before clogging occurs, clean regularly (once a week or so) or when drainage of condensed water accumulated on the door rail becomes poor. Wipe with a soft cloth when cleaning. Do not clean with a hard-bristled brush or sharp object.

Use a brush with soft bristles to remove accumulated dust.

When doing so, be careful not to damage the drainage hole and prevent clogging with dirt.

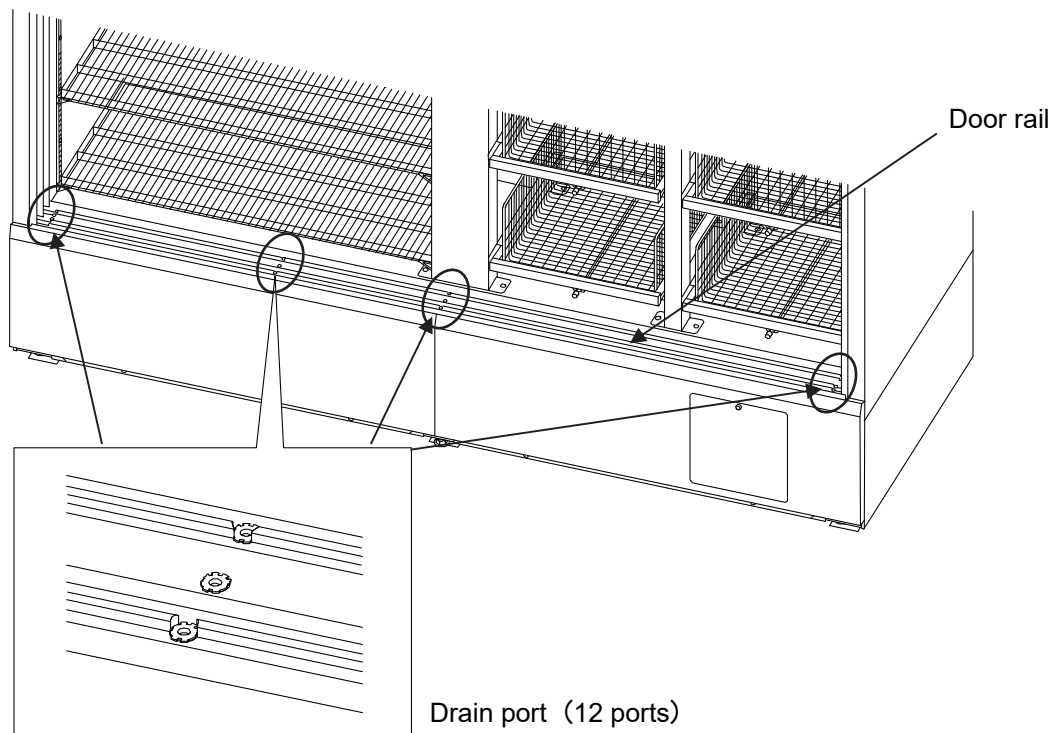

DEFROSTING

The following 2 types of defrosting methods is provided for the refrigerator. Both of them are controlled automatically.

■ Cycle defrosting

The refrigerator temperature is kept stable by ON/OFF operation of the compressor. When the compressor is OFF (stopped), the frost on the evaporator is melted by the heater. This defrosting never influences the refrigerator temperature.

■ Defrosting by defrost sensor

The cycle defrosting may not be sufficient to remove the frost on the evaporator when the ambient temperature is high, the door is opened frequently, or heavy moist items are stored in the refrigerator. In this case, the defrosting operation is started automatically when the defrost sensor detects the frost.

During the defrosting, DEF is displayed on the temperature display of home screen.

After completion of defrosting, DEF display disappears and the refrigerator returns to the normal operation.

<Important>

During the defrosting, the refrigerator temperature reaches around 10 °C temporarily.

<Reference>

Too much frost is accumulated on the evaporator when the refrigerator is running under high temperature and humidity condition. For example, the defrosting operation is started once a week when the refrigerator is running with 2 °C setting at 30 °C, 80 % R.H.

TROUBLESHOOTING

If the unit malfunctions, check the following before calling for service.

<Attention>

If the malfunction is not resolved after checking the following items or if the malfunction is not shown in the table below, contact our sales representative or agent.

Malfunction	Check/Remedy
If nothing operates even when plugged in	<ul style="list-style-type: none"> <input type="checkbox"/> The unit is not connected to the power supply properly. <input type="checkbox"/> The capacity and voltage of power supply is not sufficient. <input type="checkbox"/> There is a power failure. <input type="checkbox"/> The circuit breaker on the supply circuit is activated. <input type="checkbox"/> The fuse on the supply circuit is blown. <input type="checkbox"/> The circuit breaker switch is OFF. <input type="checkbox"/> The standby key is not pressed. <input type="checkbox"/> The plug on the unit side is not connected properly.
The alarm is activated at start-up	<ul style="list-style-type: none"> ◇ The alarm status is kept until the chamber temperature reaches the set temperature.
The alarm is activated during operation	<ul style="list-style-type: none"> <input type="checkbox"/> The unit is not connected to the power supply properly. <input type="checkbox"/> The capacity and voltage of power supply is not sufficient. <input type="checkbox"/> There is a power failure. <input type="checkbox"/> The circuit breaker on the supply circuit is activated. <input type="checkbox"/> The fuse on the supply circuit is blown. <input type="checkbox"/> The chamber temperature setting was changed. <input type="checkbox"/> The door was kept opened for a long time. <input type="checkbox"/> The items of high temperature (load) were put in the chamber. <input type="checkbox"/> The door is opened.
No key operation is available	<ul style="list-style-type: none"> <input type="checkbox"/> The key lock is set in ON. → Set the key lock in OFF [page 27].
During the setting mode, the mode returns to the temperature display mode	<ul style="list-style-type: none"> ◇ The setting mode returns to the temperature display mode automatically when 90 seconds has passed without any key operation (auto-return function).
Noisy	<ul style="list-style-type: none"> <input type="checkbox"/> The floor is not sturdy. <input type="checkbox"/> The installation site is not level. <input type="checkbox"/> The refrigerator is tilted. <input type="checkbox"/> The cabinet contact the surrounding wall.
When chamber does not get cold enough	<ul style="list-style-type: none"> <input type="checkbox"/> A large amount of items or warm product was put in the chamber. <input type="checkbox"/> The door is frequently opened. <input type="checkbox"/> The setting of chamber temperature is high. <input type="checkbox"/> The unit is in direct sunlight. <input type="checkbox"/> The unit is not installed in the installation site specified in this manual. <input type="checkbox"/> The ventilation around the unit is blocked. <input type="checkbox"/> There is a nearby heat source. <input type="checkbox"/> The ambient temperature is too high. → The allowable ambient temperature is between -5 °C and 35 °C. <input type="checkbox"/> The chamber is overloaded. <input type="checkbox"/> The air exhaust vent is blocked up with storage items. <input type="checkbox"/> The access port is not covered. → The access port should be covered with the insulation and rubber caps when no use. <input type="checkbox"/> The door gasket is damaged. → If it is damaged, contact our sales representative or agent for replacement. <input type="checkbox"/> The door cannot be fully closed..

TROUBLESHOOTING

Malfunction	Check/Remedy
Data cannot be exported to the USB flash drive (Message is displayed)	<p>“USB memory is disconnected”</p> <ul style="list-style-type: none">□ USB flash drive is not inserted into the USB port.□ USB flash drive is not inserted properly.→ Insert a USB flash drive properly into the USB port.□ USB flash drive that requires password is used.→ Use a USB flash drive that requires no password. <p>“USB memory is full”</p> <ul style="list-style-type: none">□ There is not enough free space in the USB flash drive.□ The USB flash drive is not formatted in FAT16 or FAT32 format.→ Check the capacity and format type of the USB flash drive.□ USB flash drive that requires passwords is used.→ Use a USB flash drive that requires no passwords. <p>“No Data”</p> <ul style="list-style-type: none">□ Data during the specified time period does not exist.→ Specify the time period again. <p>“An error occurred”</p> <ul style="list-style-type: none">→ Re-insert a USB flash drive or use another USB flash drive and then check the message again.□ USB flash drive that requires passwords is used.→ Use a USB flash drive that requires no passwords.

Note:

Keep an electric product which emits an electromagnetic wave away from this unit. Interference from an electromagnetic wave may cause malfunction of this unit.

DISPOSAL OF UNIT

WARNING



If the unit is to be stored unused in an unsupervised area for an extended period **ensure that children do not have access and doors cannot be closed completely.**

The disposal of the unit should be accomplished by appropriate personnel. Always remove doors to prevent accidents such as suffocation.



Flammable and explosive product. The unit contains flammable refrigerant. When repairing or recycling, only trained service personnel will repair and follow the procedure below.

- Ensure the room is well ventilated to prevent refrigerant accumulation.
- Keep fire away when the refrigerant is contained in the product.
- Do not damage or break the pipework.



Label indication is obliged to comply with Taiwanese battery regulation.

Decontamination of unit

Before disposal of unit with biological hazards, decontaminate the unit as much as possible.

TEMPERATURE RECORDER (OPTION)

The chamber temperature can be recorded and checked by installing an optional temperature recorder (MTR-G04A or MTR-0621LH).

◇Contact our sales representative or agent for the purchase of temperature recorder.


Main specifications of temperature recorder

	MTR-G04A	MTR-0621LH
Recording range	-10 °C to 40 °C	-6 °C to 20 °C
Feed speed of recorder chart	1-day/1 turn, 7-day/1 turn, 32-day/1 turn	31-day/batch
Recorder chart	Circular type	Strip type
Power source	Supplied from the refrigerator	Dry cell

◇For the installation of temperature recorder MTR-G04A, optional recorder mounting bracket MPR-S7 is necessary.

◇For the installation of temperature recorder MTR-0621LH, optional recorder mounting bracket MPR-S30W is necessary.

The figure shows the attachment location of the sensor.

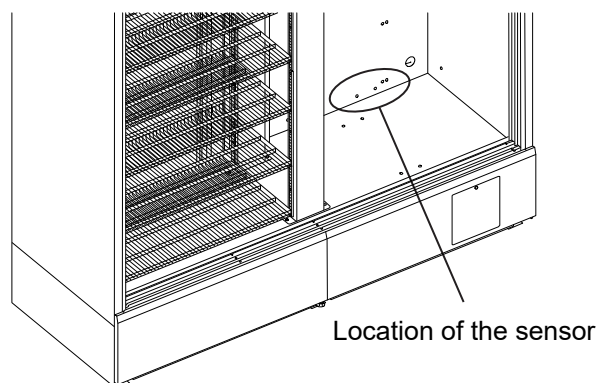

BATTERY KIT FOR POWER FAILURE ALARM (OPTION)

It is recommended to install an optional battery kit for power failure alarm in the case of power failure. At the power failure, the alarm activates to prompt an action f to prevent the stored items becoming too warm.

◇Contact our sales representative or agent for the purchase of battery kit for power failure alarm.

SPECIFICATIONS

Product name	Pharmaceutical refrigerator	
	MPR-S1201XH	MPR-S1201RXH
External dimensions	W1800 mm x D650 mm x H1824mm	
Internal dimensions	W1700 mm x D510 mm x H1425 mm	
Effective capacity	1165 L	1155 L
Exterior	Painted steel	
Interior	Painted steel	
Door	Highly insulated double glass door Glass: tempered glass	
Insulation	Rigid polyurethane foamed-in place	
Shelf	Hard steel wire with polyethylene coating x 12 Allowable load; 50 kg/shelf Inner dimensions; W781 mm x D420 mm x H23 mm	Hard steel wire with polyethylene coating x 6 Allowable load; 50 kg/shelf Inner dimensions; W781 mm x D420 mm x H23 mm
Drawer	-----	Hard steel wire powder baking finish x 10 Allowable load; 20 kg/drawer Inner dimensions; W311 mm x D355 mm x H145 mm
Access port	Inner diameter; 30 mm, 3 locations (back)	
Cooling method	Forced cool air circulation	
Compressor	Inverter type, Output; 130 W x 2	
Evaporator	Fin and tube type	
Condenser	Wire and tube type	
Refrigerant	R-600a	
Defrosting	Cycle defrost +Defrosting by defrost sensor	
Defrost heater	108 W x 2	
Drain-pan heater	25 W x 2	
Rail heater	8 W x 1	
Temperature controller	Microcomputer control system	
Temperature display	Digital display (in 1 °C, 0.1 °C increments)	
Thermal sensor	Thermistor sensor	
Memory back-up	Nonvolatile memory	
Interior light	LED x 24	
Weight	265 kg	277 kg
Accessories	2 keys, 6 side guards	2 keys, 3 side guards
Optional components	Temperature recorder (MTR-G04A, MTR-0621LH) Recorder mounting bracket (MPR-S7 for MTR-G04A) Recorder mounting bracket (MPR-S30W for MTR-0621LH) Interface board (MTR-480, MTR-L03)* Battery kit for power failure alarm (MPR-48B2) Name card holder (MPR-50CH)	

◇Design or specifications are subject to change without notice.

◇Refer to the updated catalog when ordering an optional component.

* Check the technical specifications with the sales representative or agent before purchasing.

It is recommended to use interface cables with a maximum length of 30 meters.

PERFORMANCE

Product name	Pharmaceutical Refrigerator	
	MPR-S1201XH	MPR-S1201RXH
Model number	MPR-S1201XH-PA	MPR-S1201RXH-PA
Temperature control range	2 °C to 14 °C (ambient temperature; -5 °C to 35 °C, no load)	
Noise level	42 dB (A scale)	
Maximum pressure	900 kPa	
Maximum heat emission	1404 kJ/h	
Rated voltage	AC 115 V	
Rated frequency	60 Hz	
Amperage	Max. 5.5 A	
Rated power consumption	Max. 390 W	
Usable environment condition	Temperature; -5 °C to 35 °C, Humidity; equal or less than 80 %R.H.	

◇The above data is measured based on our internal procedures.

◇Design or specifications are subject to change without notice.

SAFETY CHECK SHEET



CAUTION

Please fill in this form before servicing.
Hand over this form to the service engineer for their
and your safety.

Safety check sheet

1. Stored material

Risk of infection: Yes No Maybe

Risk of toxicity: Yes No Maybe

Risk from radioactive sources: Yes No Maybe

List all potentially hazardous materials that have been stored in this unit:

2. Contamination in the unit

a) Contamination Yes No Maybe

Types of contamination (if any): _____

b) Decontaminated Yes No

Methods used for the decontamination work:

3. Status of the unit

The unit is now safe to work on Yes No

If the answer is "No,"

Details on the danger: _____

Measures we should take to reduce the danger:

Date:

Signature:

Address, Division:

Telephone:

Product name: Pharmaceutical refrigerator	Model No. MPR	Serial number:	Date of Installation:
---	------------------	----------------	-----------------------

Please decontaminate the unit yourself before calling the service engineer.

MEMO

MEMO

MEMO



LDCL081602-1

PHC Corporation

1-1-1 Sakada, Oizumi-machi, Ora-gun, Gunma 370-0596, Japan

© PHC Corporation 2023

Verified by KABUMOTO TAKANARI 2023/10/2 16:28:50 JST

Approved by ARAI TAKASHI2 2023/10/2 16:32:28 JST

Printed in Indonesia

LDCL081602-1

N0723-10923

2023/10/3 08:23:26 JST

VALID