

Tube Rack Assembly

Model: B2000-T150/T500

The tube rack consist of 3 trays.

- The bottom tray (small holes)
- The middle tray (large holes)
- The top tray (large holes w/extra screw holes for handles).



Fig. 1

Step 1.)

Locate the two rubber coated handles and attach each to the top tray using the larger of the included screws (see Fig.2)

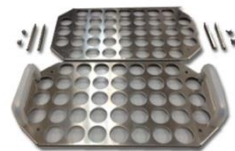


Fig. 2

Step 2)

Locate the 4 long posts & 4 small screws, attach as shown in Fig.3.



Fig. 3

Step 3)

Attach the middle tray to the other end of the long posts and thread together using the female end of the short posts. See Fig. 4 & 5



Fig. 4



Fig. 5

Step 4)

Attach the bottom tray to the short posts. See Fig. 6



Fig. 6

Step 5)

Attach the 4 plastic legs beneath the bottom plate by threading them onto the male end of the short posts. See Fig. 7



Fig. 7