

**TraceableLIVE<sup>®</sup>**  
by cole-parmer

**INTRODUCING**

# Traceable<sup>®</sup> 7600 Smart Wi-Fi Connected Data Logger

New features based on feedback from real-world labs make TraceableLIVE<sup>®</sup> more effective, more efficient, and easier to set up and use. Now you can spend more time on science and less time trying to manage your monitoring system.



**Cole-Parmer<sup>®</sup>**  
an antylia scientific company

# Traceable® 7600 Smart Wi-Fi Data Logging Refrigerator/Freezer Thermometer



## Reimagined for Today's Clinics and Labs

This new smart Wi-Fi connected data logger offers many of the best-loved features of the previous TraceableLIVE® connected data loggers, plus much more.

- Stay connected to your critical environments no matter where you are—get email, text, and push notifications to your internet-connected device
- NEW interactive and intuitive touch screen, enhanced security, and smart probes make calibration easier
- NEW acknowledge alarms remotely via the TraceableLIVE® app
- NEW calibrated smart sensor probes with easy-grip connectors allow you to quickly swap out probes for calibration and disruption-free monitoring
- NEW open-door sensors prevent accidental temperature excursions
- NEW onboard security measures offer optional PIN protection for your settings, to prevent tampering
- NEW operates using both 2.4 and 5 GHz WiFi
- Simple, one-time setup takes only a few minutes
- Includes NIST-traceable certificate of calibration from our A2LA-accredited ISO/IEC 17025:2017 (1750.01) laboratory

## Choose the Traceable® 7600 Kit That Is the Best for Your Lab

The Traceable® 7600 Smart Wi-Fi Data Logging Refrigerator/Freezer Thermometer is available in several formats so you can pick the one that fits your lab setup. The data logger and smart sensors are also available additionally. You can add to a kit or create a system for your individual needs.

## Ordering Information

Item #	Product description
07504-31	Traceable® 7600 Smart Wi-Fi Data Logging Refrigerator/Freezer Thermometer with 1 bottle probe and 1 door sensor
07504-33	Traceable® 7600 Smart Wi-Fi Data Logging Refrigerator/Freezer Thermometer with 1 bullet/ambient probe and 1 door sensor
07504-34	Traceable® 7600 Smart Wi-Fi Data Logging Refrigerator/Freezer Thermometer with 1 vaccine probe and 1 door sensor
07504-11	Traceable® 7600 Smart Wi-Fi Data Logging Refrigerator/Freezer
07504-12	Smart Sensor Calibrated Temperature Bullet Probe
07504-13	Smart Sensor Calibrated Temperature Bullet/ Ambient Probe
07504-14	Smart Sensor Calibrated Temperature Vaccine Probe
07504-29	Smart Sensor Calibrated Open-Door Sensor only

# Calibration Just Got Easier with Smart Sensors



## Keep the Data Logger and Swap the Sensor for More Efficient Recalibration

The Traceable® 7600 Smart Wi-Fi Data Logging Refrigerator/Freezer Thermometer comes equipped with smart sensor probes and open-door sensors. They're called smart because each sensor is calibrated instead of the data logger.

When it's time to recalibrate, the data logger stays put, and you can swap out the sensor that's expired for one that is ready to go. It's quick, it's easy, and it poses little to no disruption to your lab's sample monitoring activity.




## 3 Easy Steps to Recalibrate

Recalibrating is as easy as 1, 2, 3. The visual guide below will help you have your Traceable® 7600 Smart Wi-Fi Data Logging Refrigerator/Freezer Thermometer back up and running in no time.

### STEP 1

Check if calibration is near expiration.

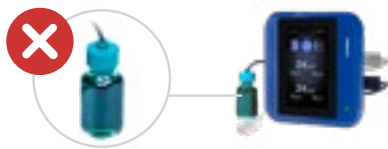

### STEP 2

Purchase new probes.



### STEP 3

Switch probes and you are ready to go.



Note: If you maintain a second set of probes, they can be returned to InnoCal for calibration.

Traceable® smart sensor probes and open-door sensors only work with the Traceable® 7600 Smart Wi-Fi Data Logging Refrigerator/Freezer Thermometer and do not operate with other data loggers. Also, noncalibrated probes or sensors are not compatible with the Traceable® 7600 Smart Wi-Fi Data Logging Refrigerator/Freezer Thermometer.

**Cole-Parmer®**  
an antylia scientific company



# TraceableLIVE® Monitors Your Samples 24/7 So You Don't Have To



## Know What's Happening in Your Refrigerator or Freezer, No Matter Where You Are

TraceableLIVE® is a cloud-based subscription software solution for connected data logging devices that sends alerts to your Wi-Fi connected device when you need to take action.

- Alerts, emails, and push notifications sent to your smartphone or Apple Watch,\* and voice call escalations\* to your phone, for peace of mind that your samples are safe
- Wi-Fi cloud-connected, so your data is available whenever you need it
- Regular updates are automatically pushed to your system, keeping it maintenance free, up to date, and secure
- Quick to set up and easy to use, keeping you focused on what matters most
- Add multiple users and multiple devices for robust lab monitoring
- Fully compliant—meets 21 CFR Part 11 North American compliance and Annex 11 for Europe and ISO 17025 accredited calibration
- The TraceableLIVE® app is kept up to date with current technological and security developments by a team of skilled programmers and professionals

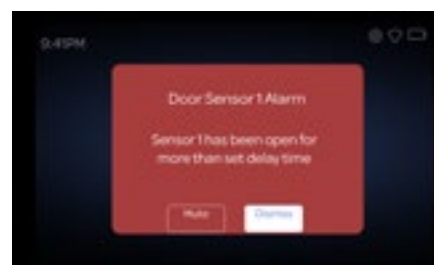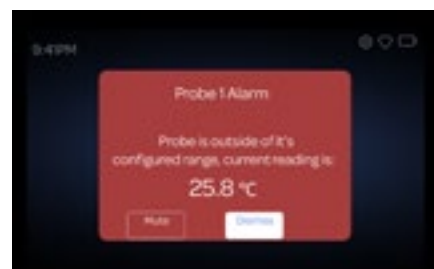

## Try TraceableLIVE® Now



Try TraceableLIVE® for yourself with our online demo or learn more at [coleparmer.com/traceablelive](https://www.coleparmer.com/traceablelive).

\*Some features and functionalities are only available with a TraceableLIVE® premium subscription. Scan the QR code to learn more.

Traceable® products have an individually numbered Traceable Certificate that assures accuracy from our A2LA-accredited ISO/IEC 17025:2017 (1750.01) calibration laboratory. It indicates traceability of measurements to the SI units through NIST or other recognized national measurement institutes (NMI) that are signatories to the CIPM Mutual Recognition Agreement. Each product goes through an intense vetting process before it can carry the Traceable name, saving you both time and money by not having to calibrate separately.

**Cole-Parmer®**  
an antylia scientific company

## TraceableLIVE® Activation and Subscription Plans

### Choose the plan to fit your facility's needs.

TraceableLIVE® is available as a monthly subscription charged to a credit card or as an annual service purchased, then initiated with an activation code.

### Subscription

Uninterrupted access to TraceableLIVE® that requires a credit card on file for monthly payments with the option of automatic annual renewal.

### Activation Code

The perfect option when using a purchase order or one-time payment that can be added to your cart and paid for a full year of TraceableLIVE® at checkout on [coleparmer.com](http://coleparmer.com). You will then receive your code and instructions via email to get started.

### TraceableLIVE® service comes in three levels.

- Free Account – Try before you buy and enjoy 30 days of Premium TraceableLIVE benefits for free.
- Standard Account – A great option for smaller teams with only a few refrigeration units to monitor.
- Premium Account – Includes helpful benefits for teams and facilities of all sizes, including escalations, recurring reports, and templates.

## Account Information

	Free Account	Standard Account	Premium Account
<b>Pricing</b>	Free	Subscription – \$8.99 per device, per month Activation – \$100 per device, per year	Subscription – \$14.99 per device, per month Activation – \$168 per device, per year
<b>Users</b>	Only one admin	Unlimited admins and users	Unlimited admins and users
<b>Max Devices</b>	3	No limit (each device must have subscription)	No limit (each device must have subscription)
<b>Data Storage</b>	Rolling 30 days	Unlimited	Unlimited
<b>Email Notifications</b>	✓	✓	✓
<b>Push-to-App Notifications</b>	✓	✓	✓
<b>Text Notifications</b>	–	–	✓
<b>Escalation Voice Call Notifications</b>	–	–	✓

Traceable® products have an individually numbered Traceable Certificate that assures accuracy from our A2LA-accredited ISO/IEC 17025:2017 (1750.01) calibration laboratory. It indicates traceability of measurements to the SI units through NIST or other recognized national measurement institutes (NMI) that are signatories to the CIPM Mutual Recognition Agreement. Each product goes through an intense vetting process before it can carry the Traceable name, saving you both time and money by not having to calibrate separately.

USA +1.800.323.4340  
+1.847.549.7600  
Canada +1.800.363.5900  
China 86.21.5109.9909  
France +33 (0) 1486 37800

Germany +49 (0) 9377 92030  
India +1.800.266.1244  
Italy +39.02.84349215  
UK +44 (0) 1480.272279  
All others +1.847.549.7600

**Cole-Parmer®**  
an antylia scientific company